




TG50R1B1 - May 10, 2022

Item # TG50R1B1 was discontinued on May 10th, 2022. For informational purposes, this is a copy of the website content at that time and is valid only for the stated product.

1X2 GRADED-INDEX (GRIN) MULTIMODE, DUAL WINDOW FIBER COUPLERS, Ø50 MM CORE

- Ø50 µm Core Graded-Index Fibers
- Dual Window Couplers: 850 nm and 1310 nm
- Available with 50:50, 75:25, 90:10, or 99:1 Coupling Ratio
- Underterminated or 2.0 mm Narrow Key FC/PC Versions

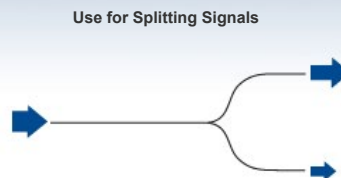

OVERVIEW

Features

- Multimode, Graded-Index (GRIN) 1x2 Couplers with Ø50 µm Core
- 850/1310 nm Dual Window Couplers
- ±40 nm Bandwidth at Specified Wavelengths
- 0.8 m Long Fiber Leads with a Tolerance of +0.075 m / -0.0 m
- Couplers Shipped with Individual Test Reports; Click Here for a Sample
- Contact Us for Custom Coupling Ratio and Connector Options



Each coupler is engraved with the item #, serial number, and key specifications for easy identification. When the white port on the left is used as the input, the coupling ratios listed below correspond to the ratio of the measured output power from the white (signal output) port to the red (tap output) port.

Thorlabs offers 1x2 graded-index (GRIN) multimode, dual-window fiber couplers with Ø50 µm core graded-index fibers. These couplers operate at both 850 ± 40 nm and 1310 ± 40 nm wavelengths. Couplers are available with 50:50, 75:25, 90:10, or 99:1 split ratios, either unterminated or with 2.0 mm narrow key FC/PC connectors. All fiber leads are 0.8 m long and are jacketed in Ø900 µm loose tubing. Additionally, each coupler ships with an individual test report; click here for a sample.

1x2 couplers are manufactured using the same process as our 2x2 fiber optic couplers, but have only one input port for simplified use and cable management. These couplers are ideal for applications where light is split from the input port into two output ports at the specified coupling ratio but are not recommended for combining applications due to losses in power. The unused port is internally terminated within the coupler housing in a manner that minimizes back reflections (please see the *1x2 Coupler Tutorial* tab for details).

These couplers are also offered with a Ø62.5 µm core size. Alternatively, Thorlabs offers standard, single-window 1x2 graded-index (GRIN) multimode fiber optic couplers with a Ø62.5 µm core.


1x2 MM Coupler Selection Guide	
Graded-Index Fiber	
Ø50 µm Core, 0.20 NA, 850 nm & 1310 nm	
Ø62.5 µm Core, 0.275 NA, 850 nm & 1310 nm	
Step-Index Fiber	
Ø50 µm Core, 0.22 NA, 400 - 900 nm (High OH)	
Ø50 µm Core, 0.22 NA, 400 - 2200 nm (Low OH)	
Ø105 µm Core, 0.22 NA, 400 - 900 nm (High OH)	
Ø105 µm Core, 0.22 NA, 400 - 2200 nm (Low OH)	
Ø200 µm Core, 0.22 NA, 400 - 900 nm (High OH)	
Ø200 µm Core, 0.22 NA, 400 - 2200 nm (Low OH)	
Ø200 µm Core, 0.39 NA, 400 - 900 nm (High OH)	
Ø200 µm Core, 0.39 NA, 400 - 2200 nm (Low OH)	
Ø200 µm Core, 0.50 NA, 400 - 900 nm (High OH)	
Ø200 µm Core, 0.50 NA, 400 - 2200 nm (Low OH)	
Ø400 µm Core, 0.39 NA, 400 - 2200 nm (Low OH)	
Ø400 µm Core, 0.50 NA, 400 - 2200 nm (Low OH)	

Alternative Fiber Coupler & Splitter Options											
Double-Clad Couplers	Single Mode Couplers			Single Mode Planar-Lightwave-Circuit Splitters		Multimode Couplers		Polarization-Maintaining Couplers			Wavelength Division Multiplexers (WDM)
	1x2	2x2	1x4	1x8	1x16	1x2	2x2	1x2	2x2	1x4	
2x2	1x2	2x2	1x4	1x8	1x16	1x2	2x2	1x2	2x2	1x4	

[Hide 1x2 Coupler Tutorial](#)

Definition of 1x2 Fused Fiber Optic Coupler Specifications

This tab provides a brief explanation of how we determine several key specifications for our 1x2 couplers. 1x2 couplers are manufactured using the same process as our 2x2 fiber optic couplers, except the second input port is internally terminated using a proprietary method that minimizes back reflections. For combining light of different wavelengths, Thorlabs offers a line of single mode wavelength division multiplexers (WDMs). The ports on our 1x2 couplers are configured as shown in the schematic below.

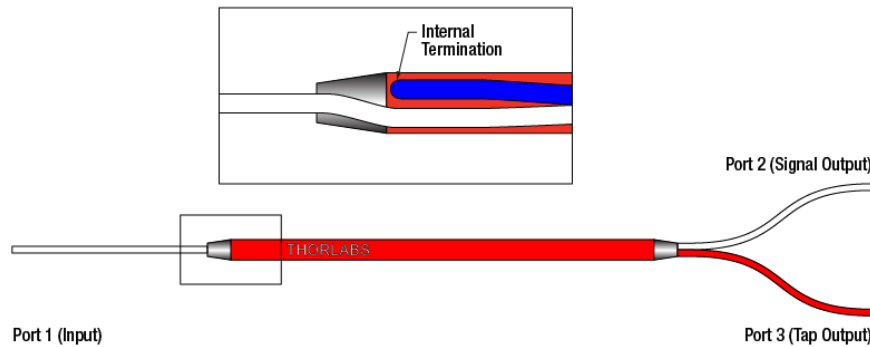

Excess Loss

Excess loss in dB is determined by the ratio of the total input power to the total output power:

$$\text{Excess Loss}(dB) = 10 \log \frac{P_{port1}(mW)}{P_{port2}(mW) + P_{port3}(mW)}$$

P_{port1} is the input power at port 1 and $P_{port2}+P_{port3}$ is the total output power from Ports 2 and 3. All powers are expressed in mW.

Polarization Dependent Loss (PDL)

The polarization dependent loss is defined as the ratio of the maximum and minimum transmissions due to polarization states in couplers. This specification pertains only to couplers not designed for maintaining polarization. PDL is always specified in decibels (dB), and can be calculated with the following equation:

$$\text{Polarization Dependent Loss}(dB) = 10 \log \frac{P_{max}(mW)}{P_{min}(mW)}$$

where P_{max} is the maximum power able to be transmitted through the coupler when scanning across all possible polarization states. P_{min} is the minimum transmission across those same states.

Optical Return Loss (ORL) / Directivity

The directivity refers to the fraction of input light that is lost in the internally terminated fiber end within the coupler housing when port 1 is used as the input. It can be calculated in units of dB using the following equation:

$$\text{Directivity}(dB) = 10 \log \frac{P_{port1b}(mW)}{P_{port1}(mW)}$$

where P_{port1} and P_{port1b} are the optical powers (in mW) in port 1 and the internally terminated fiber, respectively. This output is the result of back reflection at the junction of the legs of the coupler and represents a loss in the total light output at ports 2 and 3. For a 50:50 coupler, the directivity is equal to the optical return loss (ORL).

Insertion Loss

The insertion loss is defined as the ratio of the input power to the output power at one of the output legs of the coupler (signal or tap). Insertion loss is always specified in decibels (dB). It is generally defined using the equation below:

$$\text{Insertion Loss}(dB) = 10 \log \frac{P_{in}(mW)}{P_{out}(mW)}$$

where P_{in} and P_{out} are the input and output powers (in mW). For our 1x2 couplers, the insertion loss specification is provided for both signal and tap outputs; our specifications always list insertion loss for the signal output first. To define the insertion loss for a specific output (port 2 or port 3), the equation is rewritten as:

$$\text{Insertion Loss}_{port1 \rightarrow port2} (dB) = 10 \log \frac{P_{port1} (mW)}{P_{port2} (mW)}$$

$$\text{Insertion Loss}_{port1 \rightarrow port3} (dB) = 10 \log \frac{P_{port1} (mW)}{P_{port3} (mW)}$$

Insertion loss inherently includes both coupling (e.g., light transferred to the other output leg) and excess loss (e.g., light lost from the coupler) effects. The maximum allowed insertion loss for each output, signal and tap, are both specified. Because the insertion loss in each output is correlated to light coupled to the other output, no coupler will ever have the maximum insertion loss in both outputs simultaneously.

Calculating Insertion Loss using Power Expressed in dBm

Insertion loss can also be easily calculated with the power expressed in units of dBm. The equation below shows the relationship between power expressed in mW and dBm:


$$P(dBm) = 10 \log P(mW)$$

Then, the insertion loss in dB can be calculated as follows:

$$\text{Insertion Loss}(dB) = P_{in}(dBm) - P_{out}(dBm)$$

Coupling Ratio


Insertion loss (in dB) is the ratio of the input power to the output power from each leg of the coupler as a function of wavelength. It captures both the coupling ratio and the excess loss. The coupling ratio is calculated from the measured insertion loss. Coupling ratio (in %) is the ratio of the optical power from each output port (ports 2 and 3) to the sum of the total power of both output ports as a function of wavelength. Path A represents light traveling from port 1 to port 2 while Path B represents light traveling from port 1 to port 3. It is not impacted by spectral features such as the water absorption region because both output legs are affected equally.

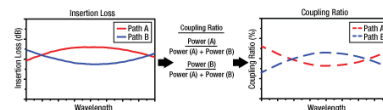

Click to Enlarge
A graphical representation of the coupling ratio calculation.

Uniformity

The uniformity is also calculated from the measured insertion loss. Uniformity is the variation (in dB) of the insertion loss over the bandwidth. It is a measure of how evenly the insertion loss is distributed over the spectral range. The uniformity of Path A is the difference between the value of highest insertion loss and the solid red insertion loss curve (in the Insertion Plot above). The uniformity of Path B is the difference between the solid blue insertion loss curve and the value of lowest insertion loss.

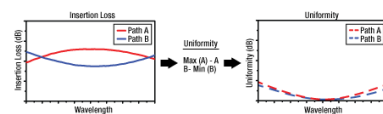

Click to Enlarge
A graphical representation of the Uniformity calculation.

[Hide Damage Threshold](#)

DAMAGE THRESHOLD


Laser-Induced Damage in Silica Optical Fibers

The following tutorial details damage mechanisms relevant to unterminated (bare) fiber, terminated optical fiber, and other fiber components from laser light sources. These mechanisms include damage that occurs at the air / glass interface (when free-space coupling or when using connectors) and in the optical fiber itself. A fiber component, such as a bare fiber, patch cable, or fused coupler, may have multiple potential avenues for damage (e.g., connectors, fiber end faces, and the device itself). The maximum power that a fiber can handle will always be limited by the lowest limit of any of these damage mechanisms.


While the damage threshold can be estimated using scaling relations and general rules, absolute damage thresholds in optical fibers are very application dependent and user specific. Users can use this guide to estimate a safe power level that minimizes the risk of damage. Following all appropriate preparation and handling guidelines, users should be able to operate a fiber component up to the specified maximum power level; if no maximum is specified for a component, users should abide by the "practical safe level" described below for safe operation of the component. Factors that can reduce power handling and cause damage to a fiber component include, but are not limited to, misalignment during fiber coupling, contamination of the fiber end face, or imperfections in the fiber itself. For further discussion about an optical fiber's power handling abilities for a specific application, please contact Thorlabs' Tech Support.

Damage at the Air / Glass Interface

There are several potential damage mechanisms that can occur at the air / glass interface. Light is incident on this interface when free-space coupling or when two fibers are mated using optical connectors. High-intensity light can damage the end face leading to reduced power handling and permanent damage to the fiber. For fibers terminated with optical connectors where the connectors are fixed to the fiber ends using epoxy, the heat generated by high-intensity light can burn the epoxy and leave residues on the fiber facet directly in the beam path.

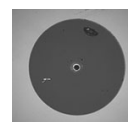

Click to Enlarge
Damaged Fiber End

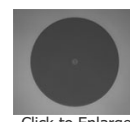

Click to Enlarge
Undamaged Fiber End

Damage Mechanisms on the Bare Fiber End Face

Estimated Optical Power Densities on Air / Glass Interface^a

Damage mechanisms on a fiber end face can be modeled similarly to bulk optics, and industry-standard damage thresholds for UV Fused Silica substrates can be applied to silica-based fiber. However, unlike bulk optics, the relevant surface areas and beam diameters involved at the air / glass interface of an optical fiber are very small, particularly for coupling into single mode (SM) fiber. therefore, for a given power density, the power incident on the fiber needs to be lower for a smaller beam diameter.

The table to the right lists two thresholds for optical power densities: a theoretical damage threshold and a "practical safe level". In general, the theoretical damage threshold represents the estimated maximum power density that can be incident on the fiber end face without risking damage with very good fiber end face and coupling conditions. The "practical safe level" power density represents minimal risk of fiber damage. Operating a fiber or component beyond the practical safe level is possible, but users must follow the appropriate handling instructions and verify performance at low powers prior to use.

Type	Theoretical Damage Threshold ^b	Practical Safe Level ^c
CW (Average Power)	~1 MW/cm ²	~250 kW/cm ²
10 ns Pulsed (Peak Power)	~5 GW/cm ²	~1 GW/cm ²

- All values are specified for unterminated (bare), undoped silica fiber and apply for free space coupling into a clean fiber end face.
- This is an estimated maximum power density that can be incident on a fiber end face without risking damage. Verification of the performance and reliability of fiber components in the system before operating at high power must be done by the user, as it is highly system dependent.
- This is the estimated safe optical power density that can be incident on a fiber end face without damaging the fiber under most operating conditions.

Calculating the Effective Area for Single Mode Fibers

The effective area for single mode (SM) fiber is defined by the mode field diameter (MFD), which is the cross-sectional area through which light propagates in the fiber; this area includes the fiber core and also a portion of the cladding. To achieve good efficiency when coupling into a single mode fiber, the diameter of the input beam must match the MFD of the fiber.

As an example, SM400 single mode fiber has a mode field diameter (MFD) of ~Ø3 µm operating at 400 nm, while the MFD for SMF-28 Ultra single mode fiber operating at 1550 nm is Ø10.5 µm. The effective area for these fibers can be calculated as follows:

$$\text{SM400 Fiber: Area} = \pi \times (\text{MFD}/2)^2 = \pi \times (1.5 \mu\text{m})^2 = 7.07 \mu\text{m}^2 = 7.07 \times 10^{-8} \text{ cm}^2$$

$$\text{SMF-28 Ultra Fiber: Area} = \pi \times (\text{MFD}/2)^2 = \pi \times (5.25 \mu\text{m})^2 = 86.6 \mu\text{m}^2 = 8.66 \times 10^{-7} \text{ cm}^2$$

To estimate the power level that a fiber facet can handle, the power density is multiplied by the effective area. Please note that this calculation assumes a uniform intensity profile, but most laser beams exhibit a Gaussian-like shape within single mode fiber, resulting in a higher power density at the center of the beam compared to the edges. Therefore, these calculations will slightly overestimate the power corresponding to the damage threshold or the practical safe level. Using the estimated power densities assuming a CW light source, we can determine the corresponding power levels as:

$$\text{SM400 Fiber: } 7.07 \times 10^{-8} \text{ cm}^2 \times 1 \text{ MW/cm}^2 = 7.1 \times 10^{-8} \text{ MW} = 71 \text{ mW (Theoretical Damage Threshold)}$$

$$7.07 \times 10^{-8} \text{ cm}^2 \times 250 \text{ kW/cm}^2 = 1.8 \times 10^{-5} \text{ kW} = 18 \text{ mW (Practical Safe Level)}$$

$$\text{SMF-28 Ultra Fiber: } 8.66 \times 10^{-7} \text{ cm}^2 \times 1 \text{ MW/cm}^2 = 8.7 \times 10^{-7} \text{ MW} = 870 \text{ mW (Theoretical Damage Threshold)}$$

$$8.66 \times 10^{-7} \text{ cm}^2 \times 250 \text{ kW/cm}^2 = 2.1 \times 10^{-4} \text{ kW} = 210 \text{ mW (Practical Safe Level)}$$

Effective Area of Multimode Fibers

The effective area of a multimode (MM) fiber is defined by the core diameter, which is typically far larger than the MFD of an SM fiber. For optimal coupling, Thorlabs recommends focusing a beam to a spot roughly 70 - 80% of the core diameter. The larger effective area of MM fibers lowers the power density on the fiber end face, allowing higher optical powers (typically on the order of kilowatts) to be coupled into multimode fiber without damage.

Damage Mechanisms Related to Ferrule / Connector Termination

Fibers terminated with optical connectors have additional power handling considerations. Fiber is typically terminated using epoxy to bond the fiber to a ceramic or steel ferrule. When light is coupled into the fiber through a connector, light that does not enter the core and propagate down the fiber is scattered into the outer layers of the fiber, into the ferrule, and the epoxy used to hold the fiber in the ferrule. If the light is intense enough, it can burn the epoxy, causing it to vaporize and deposit a residue on the face of the connector. This results in localized absorption sites on the fiber end face that reduce coupling efficiency and increase scattering, causing further damage.


For several reasons, epoxy-related damage is dependent on the wavelength. In general, light scatters more strongly at short wavelengths than at longer wavelengths. Misalignment when coupling is also more likely due to the small MFD of short-wavelength SM fiber that also produces more scattered light.

To minimize the risk of burning the epoxy, fiber connectors can be constructed to have an epoxy-free air gap between the optical fiber and ferrule near the fiber end face. Our high-power multimode fiber patch cables use connectors with this design feature.

Determining Power Handling with Multiple Damage Mechanisms

When fiber cables or components have multiple avenues for damage (e.g., fiber patch cables), the maximum power handling is always limited by the lowest damage threshold that is relevant to the fiber component. In general, this represents the highest input power that can be incident on the patch cable end face and not the coupled output power.

As an illustrative example, the graph to the right shows an estimate of the power handling limitations of a single mode fiber patch cable due to damage to the fiber end face and damage via an optical connector. The total input power handling of a terminated fiber at a given wavelength is limited by the lower of the two limitations at any given wavelength (indicated by the solid lines). A single mode fiber operating at around 488 nm is primarily limited by damage to the fiber end face (blue solid line), but fibers operating at 1550 nm are limited by damage to the optical connector (red solid line).

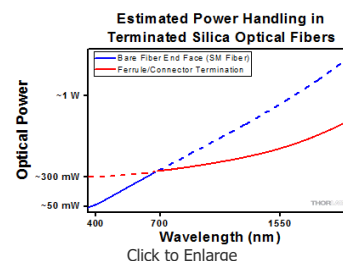

Click to Enlarge
Plot showing approximate input power that can be incident on a single mode silica optical fiber with a termination. Each line shows the estimated power level due to a specific damage mechanism. The maximum power handling is limited by the lowest power level from all relevant damage mechanisms (indicated by a solid line).

In the case of a multimode fiber, the effective mode area is defined by the core diameter, which is larger than the effective mode area for SM fiber. This results in a lower power density on the fiber end face and allows higher optical powers (on the order of kilowatts) to be coupled into the fiber without damage (not shown in graph). However, the damage limit of the ferrule / connector termination remains unchanged and as a result, the maximum power handling for a multimode fiber is limited by the ferrule and connector termination.

Please note that these are rough estimates of power levels where damage is very unlikely with proper handling and alignment procedures. It is worth noting that optical fibers are frequently used at power levels above those described here. However, these applications typically require expert users and testing at lower powers first to minimize risk of damage. Even still, optical fiber components should be considered a consumable lab supply if used at high power levels.

Intrinsic Damage Threshold

In addition to damage mechanisms at the air / glass interface, optical fibers also display power handling limitations due to damage mechanisms within the optical fiber itself. These limitations will affect all fiber components as they are intrinsic to the fiber itself. Two categories of damage within the fiber are damage from bend losses and damage from photodarkening.

Bend Losses

Bend losses occur when a fiber is bent to a point where light traveling in the core is incident on the core/cladding interface at an angle higher than the critical angle, making total internal reflection impossible. Under these circumstances, light escapes the fiber, often in a localized area. The light escaping the fiber typically has a high power density, which burns the fiber coating as well as any surrounding furcation tubing.

A special category of optical fiber, called double-clad fiber, can reduce the risk of bend-loss damage by allowing the fiber's cladding (2nd layer) to also function as a waveguide in addition to the core. By making the critical angle of the cladding/coating interface higher than the critical angle of the core/clad interface, light that escapes the core is loosely confined within the cladding. It will then leak out over a distance of centimeters or meters instead of at one localized spot within the fiber, minimizing the risk of damage. Thorlabs manufactures and sells 0.22 NA double-clad multimode fiber, which boasts very high, megawatt range power handling.

Photodarkening

A second damage mechanism, called photodarkening or solarization, can occur in fibers used with ultraviolet or short-wavelength visible light, particularly those with germanium-doped cores. Fibers used at these wavelengths will experience increased attenuation over time. The mechanism that causes photodarkening is largely unknown, but several fiber designs have been developed to mitigate it. For example, fibers with a very low hydroxyl ion (OH) content have been found to resist photodarkening and using other dopants, such as fluorine, can also reduce photodarkening.

Even with the above strategies in place, all fibers eventually experience photodarkening when used with UV or short-wavelength light, and thus, fibers used at these wavelengths should be considered consumables.

Preparation and Handling of Optical Fibers

General Cleaning and Operation Guidelines

These general cleaning and operation guidelines are recommended for all fiber optic products. Users should still follow specific guidelines for an individual product as outlined in the support documentation or manual. Damage threshold calculations only apply when all appropriate cleaning and handling procedures are followed.

1. All light sources should be turned off prior to installing or integrating optical fibers (terminated or bare). This ensures that focused beams of light are not incident on fragile parts of the connector or fiber, which can possibly cause damage.
2. The power-handling capability of an optical fiber is directly linked to the quality of the fiber/connector end face. Always inspect the fiber end prior to connecting the fiber to an optical system. The fiber end face should be clean and clear of dirt and other contaminants that can cause scattering of coupled light. Bare fiber should be cleaved prior to use and users should inspect the fiber end to ensure a good quality cleave is achieved.
3. If an optical fiber is to be spliced into the optical system, users should first verify that the splice is of good quality at a low optical power prior to high-power use. Poor splice quality may increase light scattering at the splice interface, which can be a source of fiber damage.
4. Users should use low power when aligning the system and optimizing coupling; this minimizes exposure of other parts of the fiber (other than the core) to light. Damage from scattered light can occur if a high power beam is focused on the cladding, coating, or connector.

Tips for Using Fiber at Higher Optical Power



Optical fibers and fiber components should generally be operated within safe power level limits, but under ideal conditions (very good optical alignment and very clean optical end faces), the power handling of a fiber component may be increased. Users must verify the performance and stability of a fiber component within their system prior to increasing input or output power and follow all necessary safety and operation instructions. The tips below are useful suggestions when considering increasing optical power in an optical fiber or component.

1. Splicing a fiber component into a system using a fiber splicer can increase power handling as it minimizes possibility of air/fiber interface damage. Users should follow all appropriate guidelines to prepare and make a high-quality fiber splice. Poor splices can lead to scattering or regions of highly localized heat at the splice interface that can damage the fiber.
2. After connecting the fiber or component, the system should be tested and aligned using a light source at low power. The system power can be ramped up slowly to the desired output power while periodically verifying all components are properly aligned and that coupling efficiency is not changing with respect to optical launch power.
3. Bend losses that result from sharply bending a fiber can cause light to leak from the fiber in the stressed area. When operating at high power, the localized heating that can occur when a large amount of light escapes a small localized area (the stressed region) can damage the fiber. Avoid disturbing or accidentally bending fibers during operation to minimize bend losses.
4. Users should always choose the appropriate optical fiber for a given application. For example, large-mode-area fibers are a good alternative to standard single mode fibers in high-power applications as they provide good beam quality with a larger MFD, decreasing the power density on the air/fiber interface.

5. Step-index silica single mode fibers are normally not used for ultraviolet light or high-peak-power pulsed applications due to the high spatial power densities associated with these applications.

[Hide 50:50 Fiber Optic Couplers](#)

50:50 Fiber Optic Couplers



Item # ^a	Info	Wavelength Range	Excess Loss	Insertion Loss	Directivity	NA	Core / Cladding Diameter	Fiber Type	Operating Temperature	Termination
TG50R5B1		850 ± 40 nm	<1.0 dB (Typ.)	≤4.6 dB / ≤4.6 dB	≥42.0 dB	0.20	Ø50 µm / Ø125 µm	GIF50C or Equivalent	-40 to 85 °C	No Connectors
TG50R5F1		1310 ± 40 nm								FC/PC

a. All specifications are measured without connectors during the manufacturing process.

Part Number	Description	Price	Availability
TG50R5B1	1x2 MM Dual-Window Fiber Coupler, Ø50 µm GI Fibers, 50:50 Split, No Connectors	\$110.92	Today
TG50R5F1	1x2 MM Dual-Window Fiber Coupler, Ø50 µm GI Fibers, 50:50 Split, FC/PC	\$149.73	7-10 Days

[Hide 75:25 Fiber Optic Couplers](#)

75:25 Fiber Optic Couplers



Item # ^a	Info	Wavelength Range	Excess Loss	Insertion Loss	Directivity	NA	Core / Cladding Diameter	Fiber Type	Operating Temperature	Termination
TG50R3B1		850 ± 40 nm	<1.0 dB (Typ.)	≤2.5 dB / ≤7.7 dB	≥42.0 dB	0.20	Ø50 µm / Ø125 µm	GIF50C or Equivalent	-40 to 85 °C	No Connectors
TG50R3F1		1310 ± 40 nm								FC/PC

a. All specifications are measured without connectors during the manufacturing process.

Part Number	Description	Price	Availability
TG50R3B1	1x2 MM Dual-Window Fiber Coupler, Ø50 µm GI Fibers, 75:25 Split, No Connectors	\$110.92	Today
TG50R3F1	1x2 MM Dual-Window Fiber Coupler, Ø50 µm GI Fibers, 75:25 Split, FC/PC	\$149.73	Today

[Hide 90:10 Fiber Optic Couplers](#)

90:10 Fiber Optic Couplers



Item # ^a	Info	Wavelength Range	Excess Loss	Insertion Loss	Directivity	NA	Core / Cladding Diameter	Fiber Type	Operating Temperature	Termination
TG50R2B1		850 ± 40 nm	<1.0 dB (Typ.)	≤1.6 dB / ≤11.8 dB	≥42.0 dB	0.20	Ø50 µm / Ø125 µm	GIF50C or Equivalent	-40 to 85 °C	No Connectors
TG50R2F1		1310 ± 40 nm								FC/PC

a. All specifications are measured without connectors during the manufacturing process.

Part Number	Description	Price	Availability
TG50R2B1	1x2 MM Dual-Window Fiber Coupler, Ø50 µm GI Fibers, 90:10 Split, No Connectors	\$110.92	Today
TG50R2F1	1x2 MM Dual-Window Fiber Coupler, Ø50 µm GI Fibers, 90:10 Split, FC/PC	\$149.73	Today

[Hide 99:1 Fiber Optic Couplers](#)

99:1 Fiber Optic Couplers

Item # ^a	Info	Wavelength Range	Excess Loss	Insertion Loss	Directivity	NA	Core / Cladding Diameter	Fiber Type	Operating Temperature	Termination
TG50R1B1		850 ± 40 nm	<0.5 dB (Typ.)	≤0.7 dB / ≤23.6 dB	≥42.0 dB	0.20	Ø50 µm / Ø125 µm	GIF50C or Equivalent	-40 to 85 °C	No Connectors
TG50R1F1		1310 ± 40 nm								FC/PC

a. All specifications are measured without connectors during the manufacturing process.

Part Number	Description	Price	Availability
TG50R1B1	1x2 MM Dual-Window Fiber Coupler, Ø50 µm GI Fibers, 99:1 Split, No Connectors	\$110.92	Today
TG50R1F1	1x2 MM Dual-Window Fiber Coupler, Ø50 µm GI Fibers, 99:1 Split, FC/PC	\$149.73	7-10 Days

