

WFS30-14AR - May 17, 2022

Item # WFS30-14AR was discontinued on May 17, 2022. For informational purposes, this is a copy of the website content at that time and is valid only for the stated product.

SHACK-HARTMANN WAVEFRONT SENSORS

- ▶ CMOS-Based Sensors Capable of up to 1120 fps
- ▶ Sensitivities up to $\lambda/200$
- ▶ 300 - 1100 nm or 400 - 900 nm Wavelength Range
- ▶ Kits Available with Interchangeable Microlens Arrays

Application Idea

A WFS20-5C Mounted on the KM100WFS Kinematic Mount



WFS40-7AR
 Large-Area WFS
 with 150 μ m Pitch MLA



WFS30-14AR
 General-Purpose WFS
 with 300 μ m Pitch MLA



WFS20-5C
 High-Speed WFS
 with 150 μ m Pitch MLA

OVERVIEW

Features

- Three Compact Sensor Head Options Include:
 - General-Purpose Sensor
 - Large-Aperture Sensor with 11.26 mm Square Area
 - High-Speed Sensor with Frame Rates up to 1120 fps
- High-Quality Photolithographic Microlens Array (MLA) Included
 - Mounted in Exchangeable Magnetic Holder
 - Individually Calibrated for the Sensor Head
- Kits Include Sensor Head Assembly with Additional 300 μ m Microlens Array
- Real-Time Wavefront and Intensity Distribution Measurements
- Use with CW or Pulsed Light Sources
- USB Connection for PC Operation
- Live Data Readout via TCP/IP
- Flexible Software Options
 - GUI Software
 - Instrument Driver Package for C Compilers

| Shack-Hartmann Wavefront Sensor Selection Guide ^a | | | | | |
|--------------------------------------------------------------|-----------------|------------------------|---------------------|---------------------------------|------------------------------------|
| Item # Prefix (Sensor Head) | Type | Max Speed ^b | Sensor Size | Wavefront Accuracy ^b | Wavefront Sensitivity ^b |
| WFS30 | General Purpose | Up to 64 fps | 11.34 mm x 7.13 mm | Up to $\lambda/60$ | Up to $\lambda/200$ |
| WFS40 | Large Aperture | Up to 86 fps | 11.26 mm x 11.26 mm | Up to $\lambda/40$ | Up to $\lambda/120$ |
| WFS20 | High Speed | Up to 1120 fps | 7.20 mm x 5.40 mm | Up to $\lambda/60$ | Up to $\lambda/200$ |
| Item # Suffix (MLA) ^c | Mask or Coating | Wavelength | Lenslet Pitch | Lenslet Size | Reflectivity |
| -5C | Chrome Mask | 300 - 1100 nm | 150 μ m | \varnothing 146 μ m | <25% |
| -7AR | AR Coating | 400 - 900 nm | 150 μ m | \varnothing 146 μ m | <1% |
| -14AR | AR Coating | 400 - 900 nm | 300 μ m | 295 x 295 μ m | <1% |

a. See Specs tab for complete specifications.
 b. Dependent on MLA, Operating Mode, and Specified Active Area
 c. To upgrade a wavefront sensor for compatibility with additional MLAs, contact Tech Support. Note that kits are available below which include and are calibrated for an additional -14AR MLA.

- o LabWindows™/CVI
- o LabVIEW™
- o DataSocket for Live Data Transfer
- o .Net
- Kinematic Mount Designed for Wavefront Sensors Available Separately (See Below)

Thorlabs' Fast Shack-Hartmann Wavefront Sensors provide accurate measurements of the wavefront shape and intensity distribution of incident beams. These wavefront sensors include both a CMOS-based sensor head and a microlens array (MLA) in a convenient storage and carrying case. The performance of the wavefront sensor is determined by which sensor head and MLA are combined. As summarized in the table above, we offer three sensor heads: general purpose, large aperture, and high speed. Likewise, there are three MLA options with different lenslet pitches and focal lengths for optimal accuracy and wavefront dynamic range for your application.


The Item #s for these wavefront sensors have a prefix that refers to the sensor head type (WFS30, WFS40, or WFS20), and a suffix that refers to either the MLA type (-5C, -7A, or -14AR) or the kit type (-K1 or -K2). Kits include a single sensor head and two MLAs. The *Specs* and *Selection Guide* tabs provide detailed information about the different wavefront sensor options. Trigger cables for each type of sensor head, available below, can be ordered separately for use with a TTL trigger signal. The *Shipping List* tab provides an itemized list of all components included with each item.

Sensor Heads

The general-purpose and large-aperture sensor heads, with Item # prefixes WFS30 and WFS40 respectively, have a trigger input and USB 3.0 port for connection to a PC. Operation via a USB 2.0 port will result in slower performance. High-speed sensor heads, with Item # prefix WFS20, include a control box with a trigger input and USB 2.0 port for connection to a PC. The box analyzes the camera image when the sensor is used in high-speed mode and transmits only the spot positions to the PC instead of the entire image. This lowers the volume of data sent over the USB 2.0 connection, allowing the high-speed sensors to achieve measurement speeds of >1000 fps.

Microlens Arrays

MLAs included with our wavefront sensors have a lenslet pitch of either 150 µm or 300 µm, are either chrome masked or AR Coated, and are permanently mounted in a patented (US Patent No. 8,289,504) exchangeable magnetic holder. The magnetic holder positions the microlens array to achieve correct alignment every time it is installed on the sensor head. When a kit or an additional MLA is ordered, Thorlabs includes a removal tool, which makes it easy to switch between MLAs. To purchase an additional MLA or a different combination of MLAs for use with a sensor head, please contact Tech Support. As Thorlabs must calibrate the particular sensor head for use with each specified MLA, upgrading a sensor head for use with additional MLAs requires returning the wavefront sensor to Thorlabs.

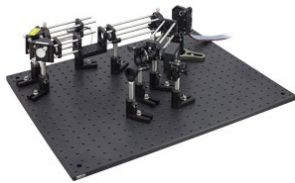

Click to Enlarge


Select wavefront sensors are available in our adaptive optics kits, shown above.

Software

The included software package offers a user-friendly graphical interface with tools for choosing camera setting, calibration, analysis, and display options. The software also includes drivers for C compilers, LabVIEW™, LabWindows/CVI™, .Net, and DataSocket server support for integration into custom system control and data collection software. For more information on the included software or to download the latest version, see the *Software* tab above.

Adaptive Optics Kits

In an effort to bring adaptive optics to even more research fields, Thorlabs provides adaptive optics kits. These kits bundle the three primary components for any adaptive optics system: a deformable mirror, a Shack-Hartmann Wavefront Sensor, and real-time control software. We offer Thorlabs' Piezoelectric Deformable Mirrors as well as MEMS-based Deformable Mirrors through our partnership with Boston Micro-Machines. In addition, the kits (shown in the photo to the left) also include a light source, all collimation/imaging optics, and all necessary mounting hardware, except for a breadboard. These kits are specifically designed to provide an affordable, easy-to-use adaptive optics solution that can be integrated into a research system in hours instead of months.

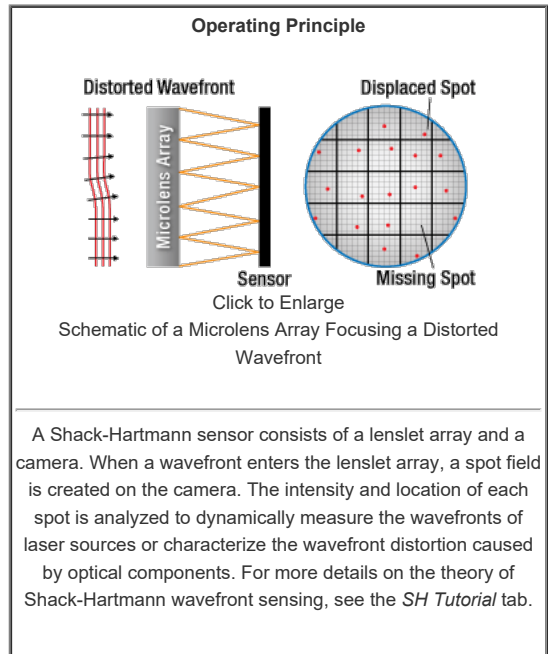

Pickup Tool for Interchanging Microlens Arrays Included with Kits

S P E C S

Thorlabs' Wavefront Sensors (WFS) combine a CMOS camera and a mounted microlens array (MLA). The performance of the wavefront sensor is primarily dictated by which WFS and MLA you choose. WFS are available with one or two MLAs. New and existing WFS can also be upgraded for use with a maximum of three types of MLA by contacting Tech Support.

The Item #s for these wavefront sensors have a prefix that refers to the sensor

| Item # Prefix | WFS30, WFS40, and WFS20 | | | | |
|-----------------------------|-------------------------|----------|-----------|---|---|
| Item # Suffix | -5C(/M) | -7AR(/M) | -14AR(/M) | - | - |
| Included Microlens Array(s) | | | | | |
| 150 µm Lenslet Pitch, | | | | | |

head type (WFS30, WFS40, or WFS20), and a suffix that refers to either the MLA type (-5C, -7A, or -14AR) or the kit type (-K1 or -K2). Specifications for the kits are not given separately. To view the specifications for a particular combination of sensor head and MLA, reference the specifications for the single-MLA Item # that includes the respective prefix and suffix.

Click the following links to move directly to the specifications of interest:

- General-Purpose WFS (Item # Prefix WFS30)
- Large-Aperture WFS (Item # Prefix WFS40)
- High-Speed WFS (Item # Prefix WFS20)

| | | | | | |
|-------------------------------------------------------------|---|---|---|---|---|
| Chrome Mask, 300 - 1100 nm | ✓ | - | - | ✓ | - |
| 150 µm Lenslet Pitch, AR ^a Coating, 400 - 900 nm | - | ✓ | - | - | ✓ |
| 300 µm Lenslet Pitch, AR ^a Coating, 400 - 900 nm | - | - | ✓ | ✓ | ✓ |

a. Anti-Reflection

General-Purpose WFS (Item # Prefix WFS30)

| Item # ^a | WFS30-5C | WFS30-7AR | WFS30-14AR |
|--------------------------------------------------------------|-------------------------------------------------------|-------------------------|--------------------------------------------|
| Microlens Arrays | | | |
| Wavelength Range | 300 nm - 1100 nm | | 400 nm - 900 nm |
| Free Aperture | 11.5 mm x 7.5 mm | | |
| Mask or Coating | Chrome Mask | Anti-Reflection Coating | |
| Reflectivity | <25% | <1% | |
| Lenslet Pitch | 150 µm | | 300 µm |
| Lenslet Grid Type | Square Grid | | |
| Lens Shape^b | Round, Plano-Convex Spherical | | Square, Plano-Convex Parabolic |
| Lens Size | Ø146 µm | | 295 µm x 295 µm |
| Fill Factor (Approximate)^c | 74.5% | | 100% |
| Number of Active Lenslets | 73 x 45 (Max), Software-Selectable Options | | 35 x 21 (Max), Software-Selectable Options |
| Effective Focal Length (Typical, When Mounted in WFS) | 4.1 mm | 5.2 mm | 14.6 mm |
| Nominal Focal Length | 5.6 mm | | 14.2 mm |
| Array Size | 12 mm x 12 mm x 1.2 mm | | |
| Substrate Material | Fused Silica (Quartz) | | |
| Camera | | | |
| Sensor Type | CMOS | | |
| Resolution | 1936 x 1216 Pixels (Max), Software-Selectable Options | | |
| Sensor Size, Max | 11.34 mm x 7.13 mm | | |
| Pixel Size | 5.83 µm x 5.83 µm | | |
| Shutter | Global ^d | | |
| Exposure Range | 33 µs - 998 ms | | |
| Image Digitization | 8 Bit | | |
| Wavefront Measurement | | | |
| Accuracy (at 633 nm)^e | λ/25 rms | λ/40 rms | λ/60 rms |
| Sensitivity (at 633 nm)^f | λ/80 rms | λ/120 rms | λ/200 rms |
| Dynamic Range (at 633 nm)^g | >500λ | >500λ | >250λ |
| Local Curvature^h | >7.4 mm | >10.0 mm | >40.0 mm |

- a. The WFS30-K1(/M) Kit operates as a WFS30-5C or WFS30-14AR, depending on which mounted MLA is installed. Similarly, the WFS20-K2(/M) Kit operates as a WFS30-7AR or WFS30-14AR.

| Software Selectable Camera Resolution Options ^a | | | | | | | |
|------------------------------------------------------------|------------------------|------------------|-----------------|------------|-------------------------------|------------------------|--|
| Camera Resolution (Pixels) | | Active Area (mm) | Microlens Spots | Spot Count | Frame Rate (fps) ^c | | |
| Normal Mode | sub2 Mode ^b | | | | Normal Mode | sub2 Mode ^b | |
| WFS30-5C, WFS30-7AR | | | | | | | |
| 1936 x 1216 | 968 x 608 | 11.34 x 7.13 | 73 x 45 | 3285 | 8.0 | 13 | |
| 1216 x 1216 | 608 x 608 | 7.13 x 7.13 | 45 x 45 | 2025 | 10 | 14 | |
| 1024 x 1024 | 512 x 512 | 6.00 x 6.00 | 39 x 39 | 1521 | 12 | 24 | |
| 768 x 768 | 384 x 384 | 4.50 x 4.50 | 29 x 29 | 841 | 21 | 32 | |
| 512 x 512 | 256 x 256 | 3.00 x 3.00 | 19 x 19 | 361 | 31 | 46 | |
| 360 x 360 | 360 x 360 | 2.11 x 2.11 | 13 x 13 | 169 | 64 | 64 | |
| WFS30-14AR | | | | | | | |
| 1936 x 1216 | 968 x 608 | 11.34 x 7.13 | 35 x 21 | 735 | 8.0 | 20 | |
| 1216 x 1216 | 608 x 608 | 7.13 x 7.13 | 21 x 21 | 441 | 13 | 20 | |
| 1024 x 1024 | 512 x 512 | 6.00 x 6.00 | 19 x 19 | 361 | 16 | 24 | |
| 768 x 768 | 384 x 384 | 4.50 x 4.50 | 13 x 13 | 169 | 21 | 32 | |
| 512 x 512 | 256 x 256 | 3.00 x 3.00 | 9 x 9 | 81 | 47 | 46 | |
| 360 x 360 | 360 x 360 | 2.11 x 2.11 | 5 x 5 | 25 | 64 | 64 | |

- a. The selected camera resolution and operation mode determines the wavefront sensor resolution (the number of calculated microlens spots) and the frame rate.
- b. When in sub2 Mode, subsampling is used to increase measurement speed. Subsampling analyzes a single pixel in each 2 x 2 block of 4 pixels.
- c. Typical speed without graphical display, depending on PC hardware, 4th order Zernike fit at specific camera resolution, and minimum exposure time.

| General Specifications | |
|----------------------------------------------|----------------------|
| External Trigger Input Specifications | |
| Trigger Slope | Software Selectable: |

- b. The convex surfaces face the camera.
- c. The Fill Factor is a measure of the fraction of incident light on the detector. For the WFS30-5C and WFS30-7AR, the fill factor is less than 100% because the round lenses are arranged on a square grid.
- d. Global shutter exposes the entire sensor at one time.
- e. Absolute accuracy using internal reference. Measured for spherical wavefronts with a known radius of curvature. Not valid for subsampling mode (sub2 Mode).
- f. Typical relative accuracy with respect to a reference wavefront (user calibration). Reference and each measurement values are averaged over 10 frames in Normal Mode.
- g. Over entire aperture of wavefront sensor.
- h. Radius of wavefront curvature over single lenslet aperture.

| | |
|----------------------------------------|--------------------------------------------------------|
| | Low-High or High-Low |
| Safe Static Input Voltage Range | 0 to 30 VDC |
| Low Level | 0.0 V to 2.0 V |
| High Level | 5.0 V to 24 V |
| Input Current (Max) | ≤10 mA |
| Pulse Width (Min) | 100 μs |
| Slew Rate (Min) | 35 V/ms |
| Housing Dimensions and Threads | |
| Optical Input Thread | Internal C-Mount |
| Dimensions (H x W x D) | 35.4 mm x 29.0 mm x 48.3 mm (1.39" x 1.14" x 1.90") |
| External Power Supply | Max 2.4 W, via USB 3.0 ^a |
| Operating Temperature | 5 to 35 °C (Non-Condensing) |
| Storage Temperature | -20 to 60 °C |
| Warm-Up Time for Rated Accuracy | 15 Minutes |
| Post Mounting^b | One 8-32 (M4) Tap, One 1/4"-20 (M6) Tap |
| Cage Mounting | - |
| Included Thread Adapter | SM1A9: External C-Mount to Internal SM1 (1.035"-40) |

- a. Operation via USB 2.0 is feasible but with reduced speed, exposure range, and power consumption.
- b. Mounting features are provided by the included removable adapter baseplate.

All technical data are valid at 23 ± 5° C and 45 ± 15% relative humidity (non-condensing).

Large-Aperture WFS (Item # Prefix WFS40)

| Item # ^a | WFS40-5C | WFS40-7AR | WFS40-14AR |
|--------------------------------------------------------------|-----------------------------------------------|-------------------------|-----------------------------------------------|
| Microlens Arrays | | | |
| Wavelength Range | 300 nm - 1100 nm | 400 nm - 900 nm | |
| Free Aperture | 11.5 mm x 11.5 mm | | |
| Mask or Coating | Chrome Mask | Anti-Reflection Coating | |
| Reflectivity | <25% | <1% | |
| Lenslet Pitch | 150 μm | 300 μm | |
| Lenslet Grid Type | Square Grid | | |
| Lens Shape^b | Round, Plano-Convex Spherical | | Square, Plano-Convex Parabolic |
| Lens Size | Ø146 μm | | 295 μm x 295 μm |
| Fill Factor (Approximate)^c | 74.5% | | 100% |
| Number of Active Lenslets | 73 x 73 (Max), Software-Selectable Options | | 35 x 35 (Max), Software-Selectable Options |
| Effective Focal Length (Typical, When Mounted in WFS) | 4.1 mm | 5.2 mm | 14.6 mm |
| Nominal Focal Length | 5.6 mm | | 14.2 mm |
| Array Size | 12 mm x 12 mm x 1.2 mm | | |
| Substrate Material | Fused Silica (Quartz) | | |
| Camera | | | |
| Sensor Type | CMOS | | |

| Software Selectable Camera Resolution Options^a | | | | | | | |
|------------------------------------------------------------------|------------------------------|-------------------------|------------------------|-------------------|-------------------------------------|------------------------------|--|
| Camera Resolution (Pixels) | | Active Area (mm) | Microlens Spots | Spot Count | Frame Rate (fps)^c | | |
| Normal Mode | sub2 Mode^b | | | | Normal Mode | sub2 Mode^b | |
| WFS40-5C, WFS40-7AR | | | | | | | |
| 2048 x 2048 | 1024 x 1024 | 11.26 x 11.26 | 73 x 73 | 5329 | 4.0 | 11 | |
| 1536 x 1536 | 768 x 768 | 8.45 x 8.45 | 55 x 55 | 3025 | 8.0 | 19 | |
| 1024 x 1024 | 512 x 512 | 5.63 x 5.63 | 35 x 35 | 1225 | 17 | 37 | |
| 768 x 768 | 384 x 384 | 4.22 x 4.22 | 27 x 27 | 729 | 26 | 55 | |
| 512 x 512 | 256 x 256 | 2.82 x 2.82 | 17 x 17 | 289 | 48 | 89 | |
| 360 x 360 | 360 x 360 | 1.98 x 1.98 | 11 x 11 | 121 | 75 | 98 | |
| WFS40-14AR | | | | | | | |
| 2048 x 2048 | 1024 x 1024 | 11.26 x 11.26 | 35 x 35 | 1225 | 5.5 | 16 | |
| 1536 x 1536 | 768 x 768 | 8.45 x 8.45 | 27 x 27 | 729 | 9.5 | 27 | |
| 1024 x 1024 | 512 x 512 | 5.63 x 5.63 | 17 x 17 | 289 | 19 | 50 | |
| 768 x 768 | 384 x 384 | 4.22 x 4.22 | 13 x 13 | 169 | 30 | 72 | |
| 512 x 512 | 256 x 256 | 2.82 x 2.82 | 7 x 7 | 49 | 57 | 98 | |
| 360 x 360 | 360 x 360 | 1.98 x 1.98 | 5 x 5 | 25 | 86 | 98 | |

- a. The selected camera resolution and operation mode determines the wavefront sensor resolution (the number of calculated microlens spots) and the frame rate.
- b. When in sub2 Mode, subsampling is used to increase measurement speed. Subsampling analyzes a single pixel in each 2 x 2 block of 4 pixels.

| | | | |
|----------------------------------------------|-------------------------------------------------------|-------------------|-------------------|
| Resolution | 2048 x 2048 Pixels (Max), Software-Selectable Options | | |
| Sensor Size, Max | 11.26 mm x 11.26 mm | | |
| Pixel Size | 5.5 μ m x 5.5 μ m | | |
| Shutter | Global ^d | | |
| Exposure Range | 77 μ s - 499 ms | | |
| Image Digitization | 8 Bit | | |
| Wavefront Measurement | | | |
| Accuracy (at 633 nm)^e | $\lambda/25$ rms | $\lambda/30$ rms | $\lambda/40$ rms |
| Sensitivity (at 633 nm)^f | $\lambda/80$ rms | $\lambda/100$ rms | $\lambda/120$ rms |
| Dynamic Range (at 633 nm)^g | >500 λ | >500 λ | >250 λ |
| Local Curvature^h | >7.4 mm | >10.0 mm | >40.0 mm |

- The WFS20-K1(/M) Kit operates as a WFS40-5C or WFS40-14AR, depending on which mounted MLA is installed. Similarly, the WFS40-K2(/M) Kit operates as a WFS40-7AR or WFS40-14AR.
- The convex surfaces face the camera.
- The Fill Factor is a measure of the fraction of incident light on the detector. For the WFS40-5C and WFS40-7AR, the fill factor is less than 100% because the round lenses are arranged on a square grid.
- Global shutter exposes the entire sensor at one time.
- Absolute accuracy using internal reference. Measured for spherical wavefronts with a known radius of curvature. Not valid for subsampling mode (sub2 Mode).
- Typical relative accuracy with respect to a reference wavefront (user calibration). Reference and each measurement values are averaged over 10 frames in Normal Mode.
- Over entire aperture of wavefront sensor.
- Radius of wavefront curvature over single lenslet aperture.

- Typical speed without graphical display, depending on PC hardware, 4th order Zernike fit at specific camera resolution, and minimum exposure time.

| General Specifications | |
|----------------------------------------------|--------------------------------------------------------|
| External Trigger Input Specifications | |
| Trigger Slope | Software Selectable: Low-High or High-Low |
| Safe Static Input Voltage Range | 0 to 30 VDC |
| Low Level | 0.0 V to 2.0 V |
| High Level | 5.0 V to 24 V |
| Input Current (Max) | \leq 10 mA |
| Pulse Width (Min) | 100 μ s |
| Slew Rate (Min) | 35 V/ms |
| Housing Dimensions and Threads | |
| Optical Input Thread | Internal C-Mount |
| Dimensions (H x W x D) | 35.4 mm x 29.0 mm x 48.3 mm (1.39" x 1.14" x 1.90") |
| External Power Supply | Max 2.9 W, via USB 3.0 ^a |
| Operating Temperature | 5 to 35 °C (Non-Condensing) |
| Storage Temperature | -20 to 60 °C |
| Warm-Up Time for Rated Accuracy | 15 Minutes |
| Post Mounting^b | One 8-32 (M4) Tap, One 1/4"-20 (M6) Tap |
| Cage Mounting | - |
| Included Thread Adapter | SM1A9: External C-Mount to Internal SM1 (1.035"-40) |

- Operation via USB 2.0 is feasible but with reduced speed, exposure range, and power consumption.
- Mounting features are provided by the included removable adapter baseplate.

All technical data are valid at 23 \pm 5° C and 45 \pm 15% relative humidity (non-condensing).

High-Speed WFS (Item # Prefix WFS20)

| Item # ^a | WFS20-5C | WFS20-7AR | WFS20-14AR |
|----------------------------------------------|-----------------------------------------------|-----------------------------------------------|------------|
| Microlens Arrays | | | |
| Wavelength Range | 300 nm - 1100 nm | 400 nm - 900 nm | |
| Free Aperture | \varnothing 9 mm | | |
| Mask or Coating | Chrome Mask | Anti-Reflection Coating | |
| Reflectivity | <25% | <1% | |
| Lenslet Pitch | 150 μ m | 300 μ m | |
| Lenslet Grid Type | Square Grid | | |
| Lens Shape^b | Round, Plano-Convex Spherical | Square, Plano-Convex Parabolic | |
| Lens Size | \varnothing 146 μ m | 295 μ m x 295 μ m | |
| Fill Factor (Approximate)^c | 74.5% | 100% | |
| Number of Active Lenslets | 47 x 35 (Max), Software-Selectable Options | 23 x 17 (Max), Software-Selectable Options | |
| Effective Focal Length | | | |

| Software Selectable Camera Resolution Options ^a | | | | | | |
|------------------------------------------------------------|------------------|-----------------|------------|-------------------------------|--------------------------------------|-----------------|
| Camera Resolution (Pixels) | Active Area (mm) | Microlens Spots | Spot Count | Frame Rate (fps) ^b | | |
| | | | | Normal Mode | Normal Mode, 2X Binning ^c | High-Speed Mode |
| WFS20-5C and WFS20-7AR | | | | | | |
| 1440 x 1080 | 7.20 x 5.40 | 47 x 35 | 1645 | 23 | 58 | 150 |
| 1080 x 1080 | 5.40 x 5.40 | 35 x 35 | 1225 | 29 | 69 | 166 |
| 768 x 768 | 3.84 x 3.84 | 23 x 23 | 529 | 50 | 144 | 350 |
| 512 x 512 | 2.56 x 2.56 | 15 x 15 | 225 | 79 | 320 | 710 |
| 360 x 360 | 1.80 x 1.80 | 11 x 11 | 121 | 116 | 630 | 880 |
| WFS20-14AR | | | | | | |
| 1440 x 1080 | 7.20 x 5.40 | 23 x 17 | 391 | 28 | 92 | 183 |
| 1080 x 1080 | 5.40 x 5.40 | 17 x 17 | 289 | 33 | 123 | 226 |
| 768 x 768 | 3.84 x 3.84 | 11 x 11 | 121 | 50 | 258 | 394 |
| 512 x 512 | 2.56 x 2.56 | 7 x 7 | 49 | 79 | 450 | 725 |
| 360 x 360 | 1.80 x 1.80 | 5 x 5 | 25 | 116 | 648 | 1120 |

- The selected camera resolution and operation mode determines

| | | | |
|----------------------------------------|-------------------------------------------------------|-------------------|-----------------|
| (Typical, When Mounted in WFS) | 4.1 mm | 5.2 mm | 14.6 mm |
| Nominal Focal Length | 5.6 mm | | 14.2 mm |
| Array Size | 10 mm x 10 mm x 1.2 mm | | |
| Substrate Material | Fused Silica (Quartz) | | |
| Camera | | | |
| Sensor Type | CMOS | | |
| Resolution | 1440 x 1080 Pixels (Max), Software-Selectable Options | | |
| Sensor Size, Max | 7.20 mm x 5.40 mm | | |
| Pixel Size | 5.0 μm x 5.0 μm | | |
| Shutter | Global ^d | | |
| Exposure Range | 4 μs - 83.3 ms | | |
| Image Digitization | 8 Bit | | |
| Wavefront Measurement | | | |
| Accuracy (at 633 nm) ^e | $\lambda/30$ rms | $\lambda/60$ rms | |
| Sensitivity (at 633 nm) ^f | $\lambda/100$ rms | $\lambda/200$ rms | |
| Dynamic Range (at 633 nm) ^g | >100 λ | | >50 λ |
| Slope (Max) ^h | $\pm 1.0^\circ$ | $\pm 0.8^\circ$ | $\pm 0.5^\circ$ |
| Local Curvature ⁱ | >7.4 mm | >10.0 mm | >40.0 mm |

- The WFS20-K1(/M) Kit operates as a WFS20-5C or WFS20-14AR, depending on which mounted MLA is installed. Similarly, the WFS20-K2(/M) Kit operates as a WFS20-7AR or WFS20-14AR.
- The convex surfaces face the camera.
- The Fill Factor is a measure of the fraction of incident light on the detector. For the WFS20-5C and WFS20-7AR, the fill factor is less than 100% because the round lenses are arranged on a square grid.
- Global shutter exposes the entire sensor at one time.
- Absolute accuracy using internal reference. Measured for spherical wavefronts with a known radius of curvature.
- Typical relative accuracy with respect to a reference wavefront (user calibration). Reference and each measurement values are averaged over 10 frames in Normal Mode.
- Over entire aperture of wavefront sensor, Normal Mode.
- Maximum local wavefront slope in High-Speed Mode.
- Radius of wavefront curvature over single lenslet aperture.

All technical data are valid at $23 \pm 5^\circ \text{C}$ and $45 \pm 15\%$ relative humidity (non-condensing).

- the wavefront sensor resolution (the number of calculated microlens spots) and the frame rate.
- Typical speed without graphical display, depending on PC hardware, 4th order Zernike fit at specific camera resolution, and minimum exposure time.
 - In Normal Mode with 2X Binning, 2 x 2 pixels are combined and averaged.

| General Specifications | |
|----------------------------------------------|--------------------------------------------------------|
| External Trigger Input Specifications | |
| Trigger Slope | Software Selectable: Low-High or High-Low |
| Safe Input Voltage Range | -0.5 to 6.5 V |
| Low Level (Max) | 1.5 V |
| High Level (Min) | 3.5 V |
| Input Impedance | >100 k Ω |
| Pulse Width (Min) | 10 μs |
| Slew Rate (Min) | 5 V/ns |
| Housing Dimensions and Threads | |
| Optical Input Thread | Internal C-Mount |
| Camera Head Dimensions (H x W x D) | 46.0 mm x 56.0 mm x 33.1 mm (1.81" x 2.20" x 1.30") |
| FPGA Box Dimensions (H x W x D) | 27.5 mm x 57.0 mm x 98.4 mm (1.08" x 2.24" x 3.87") |
| External Power Supply | 12 V, 1.5 A |
| Operating Temperature | 5 to 35 $^\circ\text{C}$ (Non-Condensing) |
| Storage Temperature | -40 to 70 $^\circ\text{C}$ |
| Warm-Up Time for Rated Accuracy | 15 Minutes |
| Post Mounting | Three 8-32 (M4) Taps |
| Cage Mounting | Four 4-40 Taps for 30 mm Cage Systems |
| Included Thread Adapter | SM1A9: External C-Mount to Internal SM1 (1.035"-40) |

SH TUTORIAL

Shack-Hartmann Wavefront Sensors

Wavefront sensors allow a user to analyze the shape of an incident beam's wavefront in order to identify or correct aberrations caused by light traveling through individual optics or optical assemblies. Shack-Hartmann wavefront sensors achieve this by dividing the beam into an array of discrete intensity points using a microlens array. These data are used to reconstruct and analyze the shape of the wavefront using Zernike polynomials. In addition to analyzing classical optics phenomena, they are increasingly used in applications where real-time monitoring of the wavefront is used to control adaptive optics with the intent of removing the wavefront distortion before creating an image.

Spot Field Overview

To measure the wavefront of a beam, the light is aligned so that it is normally incident on the microlens array at the front of the wavefront sensor. Each lenslet collects the light filling its aperture and forms a single focal spot on the CMOS camera sensor, which is located at the focal plane of the microlens array. If the wavefront is planar, all focal spots are centered directly behind each respective lens, coincident with the optical axis of each. The result is a regularly spaced grid of spots on the camera sensor, as illustrated in Figure 1. These spot locations are called reference spot positions, and they compose the reference spot field.

When the wavefront is distorted, focal spots will be displaced from their reference positions on the camera sensor. Distorted wavefronts may also result

in spot fields that include missing spots, as illustrated in Figure 2. By comparing the locations of the spots in the measured spot field with those in the reference spot field, the shape of the wavefront can be calculated.

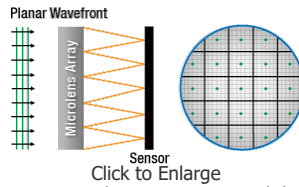


Figure 1: Focal spots are centered directly behind each lens when the incident wavefront is planar. This is the reference spot field.

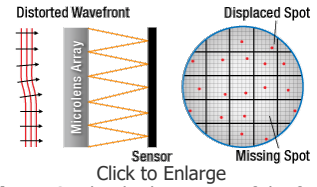


Figure 2: The displacements of the focal spots from their respective reference positions are used to calculate the shape of the incident wavefront.

Wavefront Distortion and Spot Displacement

The shape of the beam's wavefront may be complex when the entire cross section is considered; however, across the small portion of the beam collected by a single lenslet, the wavefront is approximately planar. The region surrounding a single lenslet of the microlens array is shown in Figure 3, using parallel lines to indicate both wavefronts are planar over these limited dimensions. The difference between the wavefronts is represented by their respective orientations to the optical axis of the lens.

The propagation angle of this portion of the beam is determined by the local shape of the distorted wavefront. If this region of the wavefront is not distorted, the collected light propagates along each lens' optical axis, and the focal spot is centered directly behind the lens (green circle). If instead this portion of the wavefront is distorted, its propagation angle, α , with respect to the lens' optical axis, results in a focal spot (red circle) displaced from the reference spot.

The propagation angle for this portion of the beam, and therefore the local shape of the wavefront, can be found using the focal length of the lens, f_{ML} , and the relative position of the focal spot with respect to the reference spot (δx , δy). The shape of the wavefront across the entire cross section of the beam is computed through the 2-dimensional integration of all spot displacements. Thorlabs' wavefront sensor software performs these calculations and defines the center of the spot field as the origin for the integration.

Calculating Sensitivity and Dynamic Range

Measurement Sensitivity

Sensitivity (i.e. the minimum displacement angle, α_{min}) is a function of the minimum detectable spot displacement (δy_{min}), as described by the equation:

$$\alpha_{min} = \delta y_{min} / f_{ML}$$

where f_{ML} is the focal length of the microlens.

Dynamic Range

Dynamic range (i.e. the maximum displacement angle, α_{max}) is a measure of the maximum extent of phase that can be measured:

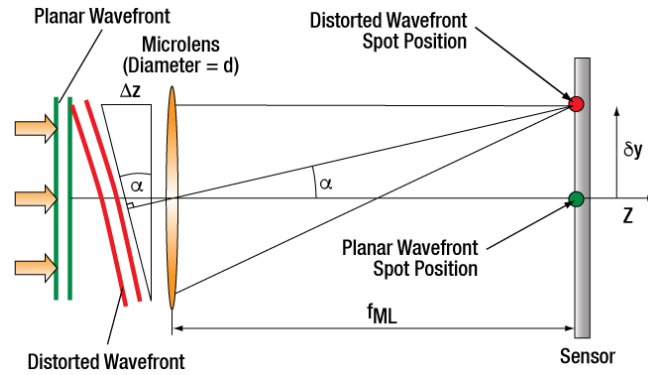


Figure 3: Above is a diagram showing a single lenslet. Waves with planar wavefronts (green parallel lines) are normally incident on the lens and focus to the reference spot position (green circle). Distorted wavefronts (red parallel lines) are not normally incident and focus to a location (red circle) displaced from the reference spot. The displacement angle, α , can be calculated from the geometry.

$$\alpha_{max} = \delta y_{max} / f_{ML} = (D / 2) / f_{ML}$$

where D is the diameter of the microlens. Both of these equations were derived using the small angle approximation. α_{min} is the minimum detectable wavefront slope that can be measured by the wavefront sensor. The minimum detectable spot displacement δy_{min} depends on the pixel size of the detector, the accuracy of the centroid algorithm, and the signal to noise ratio of the sensor. α_{max} is the maximum wavefront slope that can be measured by the wavefront sensor and corresponds to a spot displacement of δy_{max} , which is equal to the lenslet radius.

A conventional algorithm will fail to determine the correct centroid of a spot if it partially overlaps another spot or if the focal spot of a lenslet falls outside of the area of the sensor assigned to detect it (spot crossover). Special algorithms can be implemented to overcome these problems but limit the dynamic range of the sensor. The dynamic range of a system can be increased by using a lenslet with either a larger diameter or a shorter focal length. Increasing the dynamic range by increasing the lenslet diameter decreases the number of Zernike coefficients available to represent the wavefront. Conversely, increasing the dynamic range by shortening the focal length decreases the sensor's sensitivity. Ideally, a lenslet with the longest focal length that meets both the dynamic range and measurement sensitivity requirements should be used.

Factors Affecting Shack-Hartmann Measurement Accuracy

A Shack-Hartmann wavefront sensor's measurement accuracy (i.e., the minimum wavefront slope that can be measured reliably) depends on its ability to precisely measure the displacement of a focused spot with respect to a reference position. Several factors influence the measurement accuracy, including the lens pitch of the microlens array, the active area of the sensor (pupil diameter or camera resolution selection), the beam size, and the operating mode. Finer details of the wavefront can be reconstructed when there are greater numbers of focal spots to analyze. Eliminating the effects of ambient light, through a combination of software settings and experimental techniques to reduce stray light also increases measurement accuracy. Factors impacting accuracy are elaborated on below; further information can be found in the product manual.

Number of Focal Spots

The number of focal spots is determined by the pitch of the microlens array, the active area of the sensor (pupil diameter or camera resolution), and the beam size. Microlens arrays with a smaller pitch provide a greater density of measurement points for a given beam, providing higher spatial resolution, assuming all measurements are within the dynamic range of the sensor. The benefit of having larger numbers of focal spots in the spot field is particularly apparent when Zernike polynomials are used to analyze the wavefront aberrations, as the number of Zernike modes cannot exceed the number of detected focal spots. Arrays with a larger pitch provide higher wavefront accuracy.

The active area of the camera sensor is set by the user in the software. Choosing a smaller active area increases the frame rate, but information about the wavefront is lost when the active area is smaller than the beam diameter. If the beam diameter itself is small, then there will be a small number of focal spots, regardless of the active area of the camera sensor. Ideally, the beam should slightly overflow the microlens array to remove edge effects and allow any user selected area to be highlighting active pixels.

Operating Mode

Shack-Hartmann sensors can collect and average data in different ways. For our wavefront sensors, there are a total of 4 operating modes supported to varying degrees depending on the sensor type. It is important to understand that by collecting data faster, for example, may require data to be averaged. Depending on how the data is averaged, accuracy may drop. The operating modes available for our sensors are expanded on in further detail in the *Selection Guide* tab.

Image Averaging and Background Level Settings

Measurement accuracy can be increased when the average of several images, rather than a single image, is used for calculations. The wavefront sensor software includes the option to calculate normal or rolling averages. Both options compute pixel-by-pixel averages of the user-specified number of images, N. Normal averaging acquires unique sets of N images for each measurement and results in a decreased frame rate. For rolling averages, a set of N images is averaged for the first measurement, but, the next measurement will remove the first image of the set and add in the next image taken N+1. The frame rate is higher when a rolling average is used, where every frame is averaged instead of every set of frames increases. This also provides a more up-to-date status of the wavefront.

In addition to reducing stray light using lens tubes and other experimental techniques, we recommend adjusting the Noise Cut Level and Black Level settings as appropriate for the application. The software provides an intuitive and simple tool to assist with this. The Black Level setting adjusts the brightness level offset applied to all pixels. The Noise Cut Level defines a minimum brightness threshold the software uses when performing calculations, and it is used to prevent errors in focal spot calculations resulting from the effects of ambient light and/or noise. Intensity measurements below the Noise Cut Level are set to zero. The recommended Noise Cut Level setting is 'Auto,' as this setting dynamically adapts to changing light levels.

SELECTION GUIDE

Shack-Hartmann Wavefront Sensors include a sensor head and a microlens array (MLA), and the performance of the sensor depends on both. In the following, the Item # prefix of each wavefront sensor (WFS20, WFS30, or WFS40) is used to refer to the respective sensor head, while the Item # Suffixes -5C, -7AR, and -14AR are used to refer to the different MLA options. Please note that while Thorlabs offers a selection of wavefront sensors with one or two MLAs, available below, any of these sensor heads can be factory-calibrated for use with a maximum of three types of MLA by contacting Tech Support. Thorlabs also offers an upgrade service for existing WFS by contacting Tech Support.

Sensor Head

Each sensor head (Item # Prefix: WFS30, WFS40, or WFS20) is based on a different CMOS camera. While the cameras provide different measurement performance, each wavefront sensor is operated using the same software application.

Camera Responsivity

The responsivities of the three cameras, which are similar, are plotted in Figure 1. The WFS20 sensor heads have higher relative responsivities at longer wavelengths and a flatter region of peak responsivity. The WFS30 sensor has higher relative responsivities at lower wavelengths.

Free Aperture and Camera Resolution

The maximum free aperture and camera resolution are coupled for these wavefront sensors. The WFS20 and WFS30 offer rectangular apertures, with the smaller dimension limiting the maximum diameter of incident circular beams. The WFS40 is approximately 30% larger along its

| Sensor Head Selection Guide ^a | | | | | |
|------------------------------------------|-----------------|------------------------------------------|------------------------|---------------------------------|------------------------------------|
| Item # Prefix | Type | Max Sensor Size (Camera Resolution) | Max Speed ^b | Wavefront Accuracy ^b | Wavefront Sensitivity ^b |
| WFS30 | General Purpose | 11.34 mm x 7.13 mm (1936 x 1216 pixels) | Up to 64 fps | Up to $\lambda/60$ | Up to $\lambda/200$ |
| WFS40 | Large Aperture | 11.26 mm x 11.26 mm (2048 x 2048 Pixels) | Up to 289 fps | Up to $\lambda/40$ | Up to $\lambda/120$ |
| WFS20 | High Speed | 7.20 mm x 5.40 mm (1440 x 1080 pixels) | Up to 1120 fps | Up to $\lambda/60$ | Up to $\lambda/200$ |

- See Specs tab for complete specifications.
- Dependent on MLA, Operating Mode, and Specified Active Area

sides compared to the WFS20 and WFS30; it provides a square aperture that accepts an input beam of up to Ø11 mm.

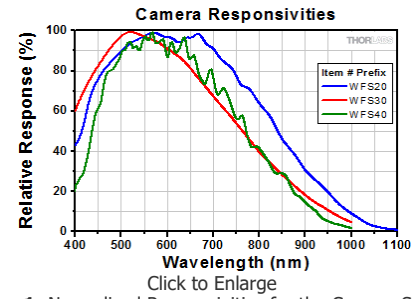


Figure 1: Normalized Responsivities for the Camera Sensors in Each Wavefront Sensor Head

Operating Modes and Measurement Speed

Wavefront sensors that include the WFS30 or WFS40 sensor heads can operate in Normal Mode or the sub2 (subsampling) Mode. There is no High-Speed Mode option offered for these sensor heads. The entire spot field image acquired by the camera is transmitted to the PC in both operating modes. When operating in sub2 Mode, the frame rate is increased by using a single pixel from each 2 x 2 set of pixels for the calculation of focal spot centroids. The drawback is wavefront measurements with reduced accuracy.

The camera included with the WFS20 operates at faster speeds than the cameras are included in the WFS30 and WFS40. Thorlabs takes advantage of the WFS20's faster camera by offering three operating modes: Normal Mode, 2 x 2 Binning Mode (Bin 2), and High-Speed Mode.

In the WFS20's normal mode, the frame rates are limited by the time needed to transmit the full spot field image from the camera to the PC, as well as by the time required to perform the calculations.

The faster 2 x 2 Binning mode improves the WFS20's frame rates by reducing calculation time. Instead of using the data from every pixel in the active area to calculate the focal spot centroids, the average value for the four pixels in every 2 x 2 set is used. The drawback is reduced spatial wavefront resolution.

In the WFS20's High-Speed Mode, the focal spot centroids are calculated by the fast field-programmable gate array (FPGA) built into the WFS20's Control Box. Not only can the FPGA calculate the spot centroids more quickly than the PC, but only the spot centroid data points are transmitted to the PC. This greatly decreases the amount of data that needs to be transferred to the PC, providing the fastest measurement speeds. A drawback is that it is not possible to view or manipulate the camera image in the GUI during operation in High-Speed Mode, which makes it harder to detect certain errors, such as those caused by camera saturation.

Physical Dimensions and Mounting Options

WFS20 sensor heads include both a sensor head and a control box, while the WFS30 and WFS40 options have only a sensor head. In addition, the dimensions of the WFS20 sensor head (46.0 mm x 56 mm x 33.1 mm) are larger than the common dimensions of the WFS30 and WFS40 sensor heads (35.4 mm x 29.0 mm x 42.6 mm).

All include an SM1A9 adapter for converting the external C-Mount to Internal SM1 (1.035"-40) threads, so that lens tubes can be used to reduce scattered light and neutral density filters can be used to prevent device saturation. The WFS20 also has three 8-32 (M4) taps on the bottom of the housing for mounting on Ø1/2" posts and four 4-40 taps on the front face for compatibility with Thorlabs' 30 mm Cage Systems. The WFS30 and WFS40 include a bottom-mounted adapter plate with an 8-32 (M4) and a 1/4"-20 (M6) tap for mounting the sensor head on Ø1/2" posts.

Microlens Arrays

The choice of microlens array affects frame rates and the quality of the wavefront measurement.

| Microlens Array Key Specifications ^a | | | | | | | | | |
|-------------------------------------------------|-----------------|---------------|---------------|-----------------|--------------|--------------------|-------------------------|-------------------------|---------------------|
| Item # Suffix | Mask or Coating | Wavelength | Lenslet Pitch | Lenslet Size | Reflectivity | High Spot Contrast | High Wavefront Accuracy | High Spatial Resolution | Low Back Reflection |
| -5C | Chrome Mask | 300 - 1100 nm | 150 µm | Ø146 µm | <25% | ✓ | - | ✓ | - |
| -7AR | AR Coating | 400 - 900 nm | 150 µm | Ø146 µm | <1% | - | - | ✓ | ✓ |
| -14AR | AR Coating | 400 - 900 nm | 300 µm | 295 µm x 295 µm | <1% | - | ✓ | - | ✓ |

a. See Specs tab for complete specifications.

Chrome Mask vs. Anti-Reflection (AR) Coating

The chrome mask is a grid that frames each microlens. The mask blocks light from being transmitted unless it passes through a microlens, which increases image contrast. However, the reflections from MLA back to the source are high. The specified wavelength range is equal to that of the quartz substrate.

The AR Coating reduces the back reflections from the MLA within the specified wavelength range. The specified reflectivity of the AR coated MLA is valid only within the stated wavelength range.


Outside of this wavelength range, the reflectivity may increase significantly. Note that the total reflectivity of the instrument may be higher due to reflections from the CMOS camera chip and its window.

Lens Pitch

An MLAs lens pitch defines the distance measured from the center of a lens to the center of the lens next to it. So, MLAs with a 150 μm lens pitch offer a larger number of spots and thus a higher spatial resolution of the wavefront, and a wider wavefront dynamic range because of their shorter focal length. The MLA with a 300 μm supports higher wavefront accuracy and sensitivity at the expense of dynamic range and spatial resolution.


Frame Rates

Frame rates depend on the sensor head, operation mode, and microlens array and scale with the number of focal spots, as is discussed in the preceding. The frame rates for each combination of sensor head, microlens array, operation mode, and spot count are plotted in Figures 2 through 4.

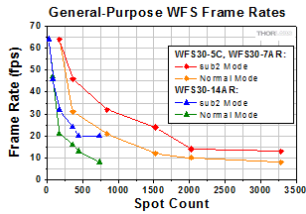


[Click to Enlarge](#)

Figure 2: Frame Rates of Wavefront Sensors with Item # Prefix WFS30

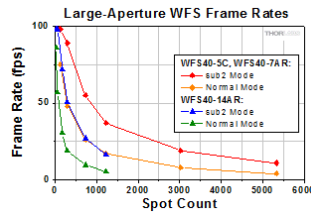

[Click to Enlarge](#)

Figure 3: Frame Rates of Wavefront Sensors with Item # Prefix WFS40

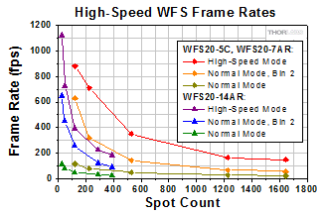

[Click to Enlarge](#)

Figure 4: Frame Rates of Wavefront Sensors with Item # Prefix WFS20

SHIPPING LIST

General-Purpose Wavefront Sensor



[Click to Enlarge](#)
WFS30-5C Package Contents

- General-Purpose Sensor Head
- Item # Specific MLA Installed on Sensor Head
 - WFS30-5C(/M): MLA with Chrome Mask and 150 μm Lenslet Pitch
 - WFS30-7AR(/M): MLA with Anti-Reflection Coating and 150 μm Lenslet Pitch
 - WFS30-14AR(/M): MLA with Anti-Reflection Coating and 300 μm Lenslet Pitch
- SM1A9 Adapter
- Dust Cap
- USB 3.0 Type-A to Mini-B Cable, 3 m
- Quick-Start Guide (Not Shown)

General-Purpose Wavefront Sensor Kit



Click to Enlarge
WFS30-K2 Package Contents

- General-Purpose Sensor Head
- Item # Specific MLA Installed on Sensor Head
 - WFS30-K1(/M): MLA with Chrome Mask and 150 μm Lenslet Pitch
 - WFS30-K2(/M): MLA with Anti-Reflection Coating and 150 μm Lenslet Pitch
- SM1A9 Adapter
- Dust Cap
- Tool to Facilitate MLA Exchange
- Additional MLA, with Anti-Reflection Coating and 300 μm Lenslet Pitch, Packaged Separately
- USB 3.0 Type-A to Mini-B Cable, 3 m
- Quick-Start Guide (Not Shown)

Large-Aperture Wavefront Sensor



Click to Enlarge
WFS40-14AR Package Contents

- Large-Aperture Sensor Head
- Item # Specific MLA Installed on Sensor Head
 - WFS40-5C(/M): MLA with Chrome Mask and 150 μm Lenslet Pitch
 - WFS40-7AR(/M): MLA with Anti-Reflection Coating and 150 μm Lenslet Pitch
 - WFS40-14AR(/M): MLA with Anti-Reflection Coating and 300 μm Lenslet Pitch
- SM1A9 Adapter
- Dust Cap
- USB 3.0 Type-A to Mini-B Cable, 3 m
- Quick-Start Guide (Not Shown)

Large-Aperture Wavefront Sensor Kit



Click to Enlarge
WFS40-K1 Package Contents

- Large-Area Sensor Head
- Item # Specific MLA Installed on Sensor Head
 - WFS40-K1(/M): MLA with Chrome Mask and 150 μm Lenslet Pitch
 - WFS40-K2(/M): MLA with Anti-Reflection Coating and 150 μm Lenslet Pitch
- SM1A9 Adapter
- Dust Cap
- Tool to Facilitate MLA Exchange
- Additional MLA, with Anti-Reflection Coating and 300 μm Lenslet Pitch, Packaged Separately
- USB 3.0 Type-A to Mini-B Cable, 3 m
- Quick-Start Guide (Not Shown)

High-Speed Wavefront Sensor



Click to Enlarge
WFS20-7AR Package Contents

- High-Speed Sensor Head
- Item # Specific MLA Installed on Sensor Head
 - WFS20-5C(/M): MLA with Chrome Mask and 150 μm Lenslet Pitch
 - WFS20-7AR(/M): MLA with Anti-Reflection Coating and 150 μm Lenslet Pitch
 - WFS20-14AR(/M): MLA with Anti-Reflection Coating and 300 μm Lenslet Pitch
- Control Box
- PoCL Cable to Connect Sensor Head and Control Box
- SM1A9 Adapter
- Dust Cap
- Two Clamps to Secure Control Box
- Power Supply with Location-Specific Plug, 12 V and 1.5 A
- USB 2.0 Type-A to Mini-B Cable, 2 m
- Quick-Start Guide (Not Shown)

High-Speed Wavefront Sensor Kit



Click to Enlarge
WFS20-7AR Package Contents

- High-Speed Sensor Head
- Item # Specific MLA Installed on Sensor Head
 - WFS20-K1(/M): MLA with Chrome Mask and 150 μm Lenslet Pitch
 - WFS20-K2(/M): MLA with Anti-Reflection Coating and 150 μm Lenslet Pitch
- Control Box
- PoCL Cable to Connect Sensor Head and Control Box
- SM1A9 Adapter
- Dust Cap
- Tool to Facilitate MLA Exchange
- Additional MLA, with Anti-Reflection Coating and 300 μm Lenslet Pitch, Packaged Separately
- Two Clamps to Secure Control Box
- Power Supply with Location-Specific Plug, 12 V and 1.5 A
- USB 2.0 Type-A to Mini-B Cable, 2 m
- Quick-Start Guide (Not Shown)

SOFTWARE

Software Download

Click on the Software button to the right for the latest version of Thorlabs' Shack-Hartmann Wavefront Sensor Software Package. The download includes the software package with a graphical user interface for operating the WFS in standard applications and support for developers who want to extend or adapt the functionality of the device to their special requirements.

System Requirements:

Software

Version 5.5 (April 9, 2021)

The wavefront sensor software package for Windows[®] 7, 8.1, 10, or later (32 or 64 bit).

- Windows® 7, 8.1, or 10 (32 or 64 Bit)
- USB Interface
 - Item # Suffix WFS20: USB 2.0 or 3.0 Port
 - Item # Suffix WFS30 or WFS40: USB 3.0 Port
- Graphic Card: 256 MB Memory (Min, Shared)
- Graphic Resolution: 1024 x 768 Min, 1280 x 1024 Recommended



Software and Graphical User Interface


Display/Output Options

For screen images of the GUI display options, please click on the links:

- Raw Spotfield Image
- Zernike Coefficients
- Measured and Reconstructed 3D Wavefront
- Irradiance Distribution
- Lineview Intensity of Pixel Column
- Tabulated Output
- Flexible Export Option: Text or Excel File
- Live Data Readout via TCP/IP to a DataSocket Server

Calculated Parameters

- Beam Centroid and Diameter
- Modal and Zonal Reconstructed Wavefront
- Max Variance of Wavefront, Peak-to-Valley (PV), and RMS of Wavefront
- Zernike Representations of Tilt, Defocus, Astigmatism, Coma, Spherical, and Higher Order Aberrations
- Fourier and Optometric Parameters

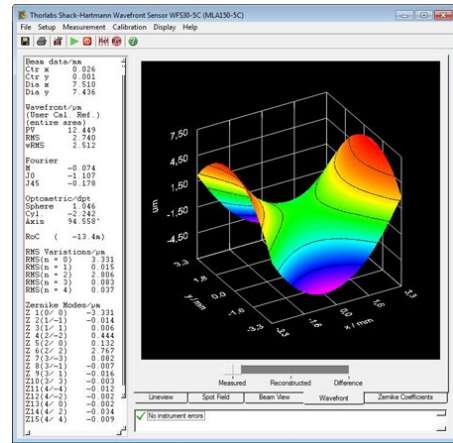

GUI Display of Measured Wavefront

Included Drivers


The software includes a driver package for constructing custom applications with the following software packages:

- C Compilers
- LabWindows™/CVI
- LabVIEW™
- .Net

General-Purpose Wavefront Sensors



- ▶ Rectangular 11.34 mm x 7.13 mm Max Aperture
- ▶ 1936 x 1216 Pixel Max Camera Resolution
- ▶ One Mounted Microlens Array Included with Each WFS
 - ▶ Chrome Mask or AR Coating
 - ▶ 150 μm Pitch Arrays Have 73 x 45 Active Lenslets (Max)
 - ▶ 300 μm Pitch Array Has 35 x 21 Active Lenslets (Max)
- ▶ User-Selectable Software Settings Control Frame Rate and Resolution
 - ▶ Pupil Size Adjustment
 - ▶ Operation in Normal or Faster Subsampling (sub2) Mode
- ▶ Software Controlled via Thorlabs' or User-Developed Application

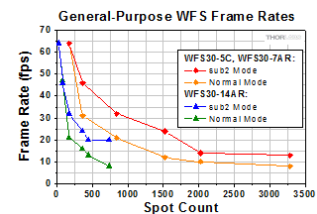

Click to Enlarge
Frame rate and resolution depend on the MLA, software-selectable active area, and operation mode.

Each of these Wavefront Sensors (WFS) is calibrated for use with the included mounted microlens array (MLA). A USB 3.0 interface to a PC running the software application is required to provide specified performance (see the Specs tab); however, these instruments can be operated with reduced performance via a USB 2.0 interface. Each WFS includes a bottom-mounted adapter plate with an 8-32 (M4) and a 1/4"-20 (M6) tap for mounting the sensor head on Ø1/2" posts.

Kits, available below, include two MLAs calibrated for the sensor head. If a MLA has not been factory-calibrated for use with a particular sensor head, it should not be used. Contact Tech Support to upgrade a WFS for use with additional MLAs.

| General Purpose Wavefront Sensors ^a | | | |
|------------------------------------------------|---------------|---------------|----------------|
| Item # | WFS30-5C(/M) | WFS30-7AR(/M) | WFS30-14AR(/M) |
| Wavelength Range | 300 - 1100 nm | | 400 - 900 nm |
| Effective Focal Length | 4.1 mm | 5.2 mm | 14.6 mm |

| | | | |
|---------------------------------------------------|---------------------------------------------|-------------------|-------------------|
| Lenslet Array Coating | Chrome Mask | AR Coated | |
| Lenslet Pitch | 150 μm | | 300 μm |
| Wavefront Accuracy (at 633 nm) | $\lambda/25$ rms | $\lambda/40$ rms | $\lambda/60$ rms |
| Wavefront Sensitivity (at 633 nm) | $\lambda/80$ rms | $\lambda/120$ rms | $\lambda/200$ rms |
| Frame Rate | 8.0 - 64 Hz | | |
| Camera Resolution | 1936 x 1216 Pixels Max, Software Selectable | | |
| High-Speed Control Box with PoCL and Table Clamps | - | | |
| Included Cable | USB 3.0 Type-A to Mini-B Cable, 3 m | | |

a. Complete specifications are available on the *Specs* tab.

| Part Number | Description | Price | Availability |
|--------------|---------------------------------------------------------------------------------------------------------------------------|------------|--------------|
| WFS30-5C/M | Shack-Hartmann WFS, 11.34 mm x 7.13 mm Aperture, 150 μm Pitch, Chrome Masked MLA, 300 - 1100 nm, Metric Taps | \$4,295.32 | Lead Time |
| WFS30-7AR/M | Shack-Hartmann WFS, 11.34 mm x 7.13 mm Aperture, 150 μm Pitch, AR Coated MLA, 400 - 900 nm, Metric Taps | \$4,295.32 | Lead Time |
| WFS30-14AR/M | Shack-Hartmann WFS, 11.34 mm x 7.13 mm Aperture, 300 μm Pitch, AR Coated MLA, 400 - 900 nm, Metric Taps | \$4,295.32 | Lead Time |
| WFS30-5C | Shack-Hartmann WFS, 11.34 mm x 7.13 mm Aperture, 150 μm Pitch, Chrome Masked MLA, 300 - 1100 nm, Imperial Taps | \$4,295.32 | Lead Time |
| WFS30-7AR | Shack-Hartmann WFS, 11.34 mm x 7.13 mm Aperture, 150 μm Pitch, AR Coated MLA, 400 - 900 nm, Imperial Taps | \$4,295.32 | Lead Time |
| WFS30-14AR | Shack-Hartmann WFS, 11.34 mm x 7.13 mm Aperture, 300 μm Pitch, AR Coated MLA, 400 - 900 nm, Imperial Taps | \$4,295.32 | Lead Time |

General-Purpose Wavefront Sensor Kits



WFS30-K1

- ▶ Includes General-Purpose Wavefront Sensor and an Additional 300 μm Pitch AR-Coated MLA
- ▶ MLAs Factory Calibrated for Use with the Included Hardware
- ▶ WFS30-K1(/M) Can Be Configured as:
 - ▶ WFS30-5C(/M) by Installing the MLA with 150 μm Pitch and Chrome Mask
 - ▶ WFS30-14AR(/M) by Installing the MLA with 300 μm Pitch and AR Coating
- ▶ WFS30-K2(/M) Can Be Configured as:
 - ▶ WFS30-7AR(/M) by Installing the MLA with 150 μm Pitch and AR Coating
 - ▶ WFS30-14AR(/M) by Installing the MLA with 300 μm Pitch and AR Coating
- ▶ Each Kit Includes a Pickup Tool Used to Easily Exchange the Installed MLA

These wavefront sensor kits come with all of the same specifications and hardware as the WFS30-5C(/M) and WFS30-7AR(/M) sold above, but they include another factory-calibrated 300 μm pitch, AR-Coated microlens array. A pickup tool is included to facilitate the easy exchange of these arrays in their patented, precision magnetic holders. Information about these wavefront sensors in the *Specs* and *Selection Guide* tabs above are given by the Item # of the configuration, rather than the Item # of the kit. For example, Item # WFS30-14AR should be referenced for the case of the 300 μm pitch MLA installed on a WFS30-K1. If a MLA has not been factory-calibrated for use with a particular sensor head, it should not be used. Contact Tech Support to upgrade a WFS for use with additional MLAs.


| Part Number | Description | Price | Availability |
|-------------|------------------------------------------------------------------------------------------------------------------------------------------------|------------|--------------|
| WFS30-K1/M | WFS Kit with 11.34 mm x 7.13 mm Aperture; 150 μm Pitch, Chrome-Masked MLA and 300 μm Pitch, AR Coated MLA; Metric Taps | \$5,165.26 | Lead Time |
| WFS30-K2/M | WFS Kit with 11.34 mm x 7.13 mm Aperture; Two AR-Coated MLAs, 150 μm Pitch and 300 μm Pitch; Metric Taps | \$5,165.26 | Lead Time |
| WFS30-K1 | WFS Kit with 11.34 mm x 7.13 mm Aperture; 150 μm Pitch, Chrome-Masked MLA and 300 μm Pitch, AR-Coated MLA; Imperial Taps | \$5,165.26 | Lead Time |
| WFS30-K2 | WFS Kit with 11.34 mm x 7.13 mm Aperture; Two AR-Coated MLAs, 150 μm Pitch and 300 μm Pitch; Imperial Taps | \$5,165.26 | Lead Time |

Large-Aperture Wavefront Sensors

- ▶ Square 11.26 x 11.26 Max Aperture
- ▶ 2048 x 2048 Pixel Max Camera Resolution
- ▶ One Mounted Microlens Array Included with Each WFS



- ▶ Chrome Mask or AR Coating
- ▶ 150 μm Pitch Arrays Have 73 x 73 Active Lenslets (Max)
- ▶ 300 μm Pitch Array Has 35 x 35 Active Lenslets (Max)
- ▶ User-Selectable Software Settings Control Frame Rate and Resolution
- ▶ Pupil Size Adjustment
- ▶ Operation in Normal or Faster Subsampling (sub2) Mode
- ▶ Software Controlled via Thorlabs' or User-Developed Application

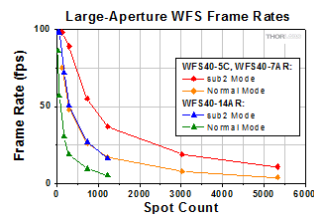

Click to Enlarge
 Frame rate and resolution (spot count) depend on the MLA, software-selectable active area, and operation mode.

Each of these Wavefront Sensors (WFS) is calibrated for use with the included mounted microlens array (MLA). A USB 3.0 interface to a PC running the software application is required to provide specified performance (see the *Specs* tab); however, these instruments can be operated with reduced performance via a USB 2.0 interface. Each WFS includes a bottom-mounted adapter plate with an 8-32 (M4) and a 1/4"-20 (M6) tap for mounting the sensor head on $\varnothing 1/2$ " posts.

Kits, available below, include two MLAs calibrated for the sensor head. If a MLA has not been factory-calibrated for use with a particular sensor head, it should not be used. Contact Tech Support to upgrade a WFS for use with additional MLAs.

| Large-Aperture Wavefront Sensors ^a | | | |
|---------------------------------------------------|---------------------------------------------|-------------------|-------------------|
| Item # | WFS40-5C(/M) | WFS40-7AR(/M) | WFS40-14AR(/M) |
| Wavelength Range | 300 - 1100 nm | 400 - 900 nm | |
| Effective Focal Length | 4.1 mm | 5.2 mm | 14.6 mm |
| Lenslet Array Coating | Chrome Mask | AR Coated | |
| Lenslet Pitch | 150 μm | | 300 μm |
| Wavefront Accuracy (at 633 nm) | $\lambda/25$ rms | $\lambda/30$ rms | $\lambda/40$ rms |
| Wavefront Sensitivity (at 633 nm) | $\lambda/80$ rms | $\lambda/100$ rms | $\lambda/120$ rms |
| Frame Rate | 4.0 - 98 Hz | | 5.5 - 98 Hz |
| Camera Resolution | 2048 x 2048 Pixels Max, Software Selectable | | |
| High-Speed Control Box with PoCL and Table Clamps | - | | |
| Included Cable | USB 3.0 Type-A to Mini-B Cable, 3 m | | |

a. Complete specifications are available on the *Specs* tab.

| Part Number | Description | Price | Availability |
|--------------|------------------------------------------------------------------------------------------------------------------------|------------|--------------|
| WFS40-5C/M | Shack-Hartmann WFS, 11.26 mm Square Aperture, 150 μm Pitch, Chrome Masked MLA, 300 - 1100 nm, Metric Taps | \$5,763.34 | Lead Time |
| WFS40-7AR/M | Shack-Hartmann WFS, 11.26 mm Square Aperture, 150 μm Pitch, AR Coated MLA, 400 - 900 nm, Metric Taps | \$5,763.34 | Lead Time |
| WFS40-14AR/M | Shack-Hartmann WFS, 11.26 mm Square Aperture, 300 μm Pitch, AR Coated MLA, 400 - 900 nm, Metric Taps | \$5,763.34 | Lead Time |
| WFS40-5C | Shack-Hartmann WFS, 11.26 mm Square Aperture, 150 μm Pitch, Chrome Masked MLA, 300 - 1100 nm, Imperial Taps | \$5,763.34 | Lead Time |
| WFS40-7AR | Shack-Hartmann WFS, 11.26 mm Square Aperture, 150 μm Pitch, AR Coated MLA, 400 - 900 nm, Imperial Taps | \$5,763.34 | Lead Time |
| WFS40-14AR | Shack-Hartmann WFS, 11.26 mm Square Aperture, 300 μm Pitch, AR Coated MLA, 400 - 900 nm, Imperial Taps | \$5,763.34 | Lead Time |

Large-Aperture Wavefront Sensor Kits



- ▶ Includes Large-Aperture Wavefront Sensor and an Additional 300 μm Pitch AR-Coated MLA
- ▶ MLAs Factory Calibrated for Use with the Included Hardware
- ▶ WFS40-K1(/M) Can Be Configured as:
 - ▶ WFS40-5C(/M) by Installing the MLA with 150 μm Pitch and Chrome Mask
 - ▶ WFS40-14AR(/M) by Installing the MLA with 300 μm Pitch and AR Coating
- ▶ WFS40-K2(/M) Can Be Configured as:
 - ▶ WFS40-7AR(/M) by Installing the MLA with 150 μm Pitch and AR Coating
 - ▶ WFS40-14AR(/M) by Installing the MLA with 300 μm Pitch and AR Coating
- ▶ Each Kit Includes a Pickup Tool Used to Easily Exchange the Installed MLA

These wavefront sensor kits come with all of the same specifications and hardware as the WFS40-5C(/M) and WFS40-7AR(/M) sold above, but they include another factory-calibrated 300 μm pitch, AR-Coated microlens array and a tool to facilitate the easy exchange of these arrays in their patented, precision magnetic holders. Information about these wavefront sensors in the *Specs* and *Selection Guide* tabs above are given by the Item # of the configuration, rather than the Item # of


the kit. For example, Item # WFS40-14AR should be referenced for the case of the 300 μm pitch MLA installed on a WFS40-K1. If a MLA has not been factory-calibrated for use with a particular sensor head, it should not be used. Contact Tech Support to upgrade a WFS for use with additional MLAs.

| Part Number | Description | Price | Availability |
|-------------|-----------------------------------------------------------------------------------------------------------------------|------------|--------------|
| WFS40-K1/M | WFS Kit with 11.26 mm Square Aperture; 150 μm Pitch, Chrome-Masked MLA and 300 μm Pitch, AR-Coated MLA; Metric Taps | \$6,633.28 | Lead Time |
| WFS40-K2/M | WFS Kit with 11.26 mm Square Aperture; Two AR-Coated MLAs, 150 μm Pitch and 300 μm Pitch; Metric Taps | \$6,633.28 | Lead Time |
| WFS40-K1 | WFS Kit with 11.26 mm Square Aperture; 150 μm Pitch, Chrome-Masked MLA and 300 μm Pitch, AR-Coated MLA; Imperial Taps | \$6,633.28 | Lead Time |
| WFS40-K2 | WFS Kit with 11.26 mm Square Aperture; Two AR-Coated MLAs, 150 μm Pitch and 300 μm Pitch; Imperial Taps | \$6,633.28 | Lead Time |

High-Speed Wavefront Sensors



- ▶ Rectangular 7.20 x 5.40 mm Max Aperture
- ▶ 1440 x 1080 Pixel Max Camera Resolution
- ▶ One Mounted Microlens Array Included with Each WFS
 - ▶ Chrome Mask or AR Coating
 - ▶ 150 μm Pitch Arrays Have 47 x 35 Active Lenslets (Max)
 - ▶ 300 μm Pitch Array Has 23 x 17 Active Lenslets (Max)
- ▶ User-Selectable Software Settings Control Frame Rate and Resolution
 - ▶ Pupil Size Adjustment
 - ▶ Operation in Normal, 2x2 Binning (Bin 2), or High-Speed Mode
- ▶ Software Controlled via Thorlabs' or User-Developed Application

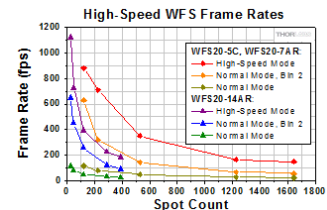

Click to Enlarge
Frame rate and resolution depend on the MLA, software-selectable active area, and operation mode.

These WFS include a separate control box that connects to the sensor head with the included cable. During high-speed mode operation, the field-programmable gate array (FPGA) built into the control box processes the image data and sends only the calculated spot centroid coordinates to the PC. In contrast, the entire image is sent to the PC when the WFS operates in normal and Bin 2 modes. A USB 2.0 interface to a PC running the software application is required for operation. Each WFS has three 8-32 (M4) taps on the bottom of the housing for mounting on Ø1/2" posts and four 4-40 taps on the front face for compatibility with Thorlabs' 30 mm Cage Systems.

Kits, available below, include two MLAs calibrated for the sensor head. If a MLA has not been factory-calibrated for use with a particular sensor head, it should not be used. Contact Tech Support to upgrade a WFS for use with additional MLAs.

| High-Speed Wavefront Sensors ^a | | | |
|---------------------------------------------------|---------------------------------------------|---------------|----------------|
| Item # | WFS20-5C(/M) | WFS20-7AR(/M) | WFS20-14AR(/M) |
| Wavelength Range | 300 - 1100 nm | 400 - 900 nm | |
| Effective Focal Length | 4.1 mm | 5.2 mm | 14.6 mm |
| Lenslet Array Coating | Chrome Mask | AR Coated | |
| Lenslet Pitch | 150 μm | | 300 μm |
| Wavefront Accuracy (at 633 nm) | λ/30 rms | | λ/60 rms |
| Wavefront Sensitivity (at 633 nm) | λ/100 rms | | λ/200 rms |
| Frame Rate | 23 - 880 Hz | | 28 - 1120 Hz |
| Camera Resolution | 1440 x 1080 Pixels Max, Software Selectable | | |
| High-Speed Control Box with PoCL and Table Clamps | ✓ | | |
| Included Cable | USB 2.0 Type-A to Mini-B Cable, 2 m | | |

a. Complete specifications are available on the [Specs](#) tab.

| Part Number | Description | Price | Availability |
|--------------|---------------------------------------------------------------------------------|------------|--------------|
| WFS20-5C/M | Fast Shack-Hartmann WFS, 150 μm Pitch, Chrome Masked, 300 - 1100 nm, M4 Taps | \$6,699.39 | Today |
| WFS20-7AR/M | Fast Shack-Hartmann WFS, 150 μm Pitch, AR Coated: 400 - 900 nm, M4 Taps | \$6,699.39 | Lead Time |
| WFS20-14AR/M | Fast Shack-Hartmann WFS, 300 μm Pitch, AR Coated: 400 - 900 nm, M4 Taps | \$6,699.39 | Today |
| WFS20-5C | Fast Shack-Hartmann, WFS, 150 μm Pitch, Chrome Masked, 300 - 1100 nm, 8-32 Taps | \$6,699.39 | Today |
| WFS20-7AR | Fast Shack-Hartmann WFS, 150 μm Pitch, AR Coated: 400 - 900 nm, 8-32 Taps | \$6,699.39 | 7-10 Days |

High-Speed Wavefront Sensor Kits



- ▶ Includes a High-Speed Wavefront Sensor, an Additional 300 µm Pitch AR-Coated MLA, and Control Box
- ▶ MLAs Factory Calibrated for Use with the Included Hardware
- ▶ WFS20-K1(/M) Can Be Configured as:
 - ▶ WFS20-5C(/M) by Installing the MLA with 150 µm Pitch and Chrome Mask
 - ▶ WFS20-14AR(/M) by Installing the MLA with 300 µm Pitch and AR Coating
- ▶ WFS20-K2(/M) Can Be Configured as:
 - ▶ WFS20-7AR(/M) by Installing the MLA with 150 µm Pitch and AR Coating
 - ▶ WFS20-14AR(/M) by Installing the MLA with 300 µm Pitch and AR Coating
- ▶ Each Kit Includes a Pickup Tool Used to Easily Exchange the Installed MLA

These wavefront sensor kits come with all of the same specifications and hardware as the WFS20-5C(/M) and WFS20-7AR(/M) sold above, but they include another factory-calibrated 300 µm pitch, AR-Coated microlens array and a tool to facilitate the easy exchange of these arrays in their patented, precision magnetic holders. Information about these wavefront sensors in the *Specs* and *Selection Guide* tabs above are given by the Item # of the configuration, rather than the Item # of the kit. For example, Item # WFS20-14AR should be referenced for the case of the 300 µm pitch MLA installed on a WFS20-K1. If a MLA has not been factory-calibrated for use with a particular sensor head, it should not be used. Contact Tech Support to upgrade a WFS for use with additional MLAs.

| Part Number | Description | Price | Availability |
|-------------|-------------------------------------------------------------------------------------------------------------|------------|--------------|
| WFS20-K1/M | Fast WFS Kit with 150 µm Pitch, Chrome-Masked and 300 µm Pitch, AR-Coated Microlens Arrays, Metric, M4 Taps | \$7,608.91 | Today |
| WFS20-K2/M | Fast WFS Kit with 150 µm Pitch and 300 µm Pitch AR-Coated Microlens Arrays, M4 Taps | \$7,608.91 | Today |
| WFS20-K1 | Fast WFS Kit with 150 µm Pitch, Chrome-Masked and 300 µm Pitch, AR-Coated Microlens Arrays, 8-32 Taps | \$7,608.91 | 7-10 Days |
| WFS20-K2 | Fast WFS Kit with 150 µm Pitch and 300 µm Pitch AR-Coated Microlens Arrays, 8-32 Taps | \$7,608.91 | Lead Time |

USB and Trigger Cables

The CAB-WFS20-T1 cable is designed to connect a user-supplied external trigger source to the High-Speed Shack-Hartmann Wavefront Sensors (Item # Prefix WFS20). The CAB-DCU-T3 cable provides the same functionality for the General-Purpose and Large-Aperture Shack-Hartmann Wavefront Sensors (Item # Prefix WFS30 and WFS40, respectively). For trigger input specifications, please see the *Specs* tab above.

| Item # | CAB-WFS20-T1 | | | Pin | Assignment |
|----------------------|-----------------------|----------------------|--|-----|-----------------------------------------|
| Click to Enlarge | Connector Device Side | 4-Pin Hirose, Female | | 1 | Not Used |
| | TTL Trigger Input | Bare Wire (Brown) | | 2 | Trigger Input Ground (Black, Bare Wire) |
| | Shielding | None | | 3 | TTL Trigger Input (Brown, Bare Wire) |
| | Length | 2 m | | 4 | Not Used |

| Item # | CAB-DCU-T3 | | | Pin | Assignment |
|----------------------|---------------------------------|----------------------------------------------------------|--|-----|--------------------------------------------|
| Click to Enlarge | Connector Device Side | Hirose HR25-7TP-8S, Female | | 1 | Ground (Gray) |
| | End Opposite Connectors | Tinned End of Wires | | 2 | Flash Output ^a - (Green) |
| | Function | GPIO | | 3 | GPIO 1, 3.3 V LVCMOS (Blue) |
| | Trigger In (Bare Wire) | Yes | | 4 | Trigger Input ^a - (Brown) |
| | Flash & Digital Out (Bare Wire) | Yes | | 5 | Flash Output ^a + (Yellow) |
| | Cable Type | Shielded High-Flexible Control Cable 8 x 0.1 mm, Ø4.9 mm | | 6 | GPIO 2, 3.3 V LVCMOS (Red) |
| | Shielding | Single Shielded | | 7 | Trigger Input ^a + (White) |
| | Length | 2 m | | 8 | Output Supply Voltage, 5 V (100 mA) (Pink) |

a. These pins are opto-decoupled inside the camera to protect against high or incorrect voltages.

| Part Number | Description | Price | Availability |
|-------------|-------------|-------|--------------|
| | | | |

| | | | |
|--------------|------------------------------------------------------------------------------|----------|-----------|
| CAB-WFS20-T1 | Trigger Cable for WFS20 Wavefront Sensors | \$78.20 | 7-10 Days |
| CAB-DCU-T3 | Trigger and I/O Cable, Hirose 25, for DCC3240, DCC3260, WFS30 and WFS40, 2 m | \$106.41 | Today |

Kinematic Mount for Wavefront Sensors



Click to Enlarge
The KM200PM mount platform's larger size allows for flexible mounting options.

- ▶ Two 80 TPI Adjusters for $\pm 4^\circ$ of Fine Tip and Tilt Control
- ▶ 8 mrad Adjustment per Revolution
- ▶ Platform can be Removed and Secured to Either Arm Enabling Left- or Right-Handed Configuration
- ▶ Removable Knobs Expose Adjuster Screws with Hex Socket

Thorlabs' KM100WFS and KM200PM(/M) Kinematic Platform Mounts provide kinematic control for our CMOS-Based Shack-Hartmann Wavefront Sensors. Please note that the optical axis of the sensors will not be centered over the optical post. The animation to the right shows

how to convert between the left- and right-handed orientations for the KM100WFS. This process is very similar for the KM200PM(/M).

The KM100WFS is designed to accommodate the WFS20 wavefront sensors, and features a similar design to our KM100PM platform. The WFS20 high-speed sensors can be mounted using the two #8 through holes near the front edge of the plate and the included 8-32 screws. The DCU CCD Cameras and the WFS20 sensors can also be mounted on the KM100WFS.

The KM200PM(/M) mount is larger than the KM100WFS, and can mount the WFS20 sensors using the two #8 (M4) through holes near the front edge of the plate. WFS30 and WFS40 can be mounted using one of the 8-32 tapped mounting holes. The larger profile and number of holes provides a variety of mounting options. The platform is secured to the front plate using two 4-40 screws with a 3/32" hex [the KM200PM/M uses M2.5 screws with a 2 mm (5/64") hex]. The two pieces of the mounting platform assembly are held together with two 3-48 screws [5/64" (2 mm) hex]. If desired, the mounting plate can also be removed and replaced with a user-supplied alternative, enabling custom mounting options.

| Platform Specifications | | | |
|-------------------------|-------------------------|--------------------|-----------------------------|
| Item # | Tapped Holes | Through Holes | Platform Depth ^a |
| KM100WFS | - | #8 (x2) M3 (x3) | 1.52" (38.6 mm) |
| KM200PM | 6-32 (x16) 8-32 (x4) | #8 (x2) | 2.10" (53.3 mm) |
| KM200PM/M | M4 (x21) | M4 (x2) | 53.3 mm (2.10") |

- a. The Distance between the Front and Back of the Platform

| Part Number | Description | Price | Availability |
|-------------|-------------------------------------------------------|----------|--------------|
| KM200PM/M | Kinematic Prism Mount, 54.4 mm Deep, M4 Taps | \$130.88 | Today |
| KM100WFS | Kinematic Mount for Thorlabs' Wavefront Sensors | \$90.68 | Today |
| KM200PM | Kinematic Prism Mount, 2.14" Deep, 6-32 and 8-32 Taps | \$130.88 | Today |

